PLT connect

Software

Impressum

BRAND GMBH + CO KG
Otto-Schott-Str. 25
97877 Wertheim (Germany)
T +49 9342 808 0
F +49 9342 808 9800
info@brand.de
www.brand.de

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Contents

1	Introduction	19
1.1	PLT unit.....	19
1.2	PLT connect.....	20
2	Install programm	20
2.1	System requirements.....	20
2.2	Rights.....	20
2.3	Before installing.....	21
2.4	Installing PLTconnect	21
3	Connecting the PLT unit to the PC.....	21
4	Operate the program.....	22
4.1	Connect to PLT unit	22
4.2	User settings.....	22
4.3	Creating pipettes.....	24
4.4	Carrying out testing	25
4.5	Filter function.....	28
4.6	Import/export data	29
4.7	Self-test	30
5	Markings	30
6	Disposal	31

1 Introduction

Before using the PLTconnect software, please read through these operating instructions as well as the instructions for the PLT unit carefully.

1.1 PLT unit

Air cushion pipettes must be checked at regular intervals within the scope of inspection equipment monitoring and the results compared with the error limits stipulated by ISO 8655-2. However, calibration certificates only provide the results from the time of the testing. The periods between these calibration procedures is critical as leaks can occur at any time.

Significantly more than 80% of pipettes returned for repair have leaks and are outside the volumetric tolerances, even although they do not drip. The most common cause for the lack of accuracy in piston-stroke pipettes is system leaks. Leaks arise through damage to the seals, pistons or tip sampling cones. Often invisible to the eye, these faults can lead to significant volumetric errors.

The PLT unit from BRAND, the leak-checking device for air cushion pipettes, detects the smallest of leaks in seconds!

1.2 PLT connect

The PLT unit software, also known as PLTconnect, is a program for facilitating the documentation of leak checks carried out with the PLT unit from BRAND. The test results are transferred from the PLT unit into the program by means of automatic measured value transmission. The result is displayed as a test certificate and saved in the program. Documentation of your test results is made significantly easier.

The program supports leak checking with and without tips as well as static and dynamic measurements.

Master data and limit values for standard single-channel and multi-channel pipettes in volumetric ranges from 1 µl to 10 ml are already stored in the program for your use. Testing is independent of the device manufacturer. You can also easily create new devices.

2 Install programm

2.1 System requirements

In order to be able to use the operating software, your PC or laptop must meet the following minimum system requirements:

- Windows® 7 32-Bit and 64-Bit
- USB 2.0 interface
- Microsoft® Installer 3.1
- NET-Framework Version 4

2.2 Rights

You require software installation rights and the necessary write rights for your PC. If you do not have these rights, contact your system administrator.


2.3 Before installing

If there is an older version of the PLT unit software already installed on your PC, this should be uninstalled first. Any log data already existing on your system will be retained after the uninstall process.

2.4 Installing PLTconnect

- a. Start "setup.exe" from the supplied USB stick and follow the setup wizard.
 - ⇒ If you accept the standard settings, the program is installed in C: \Program Files(x86)\BRAND\PLTconnect.
 - ⇒ After the installation has been successfully completed, the icon to start the program can be found on the desktop.

3 Connecting the PLT unit to the PC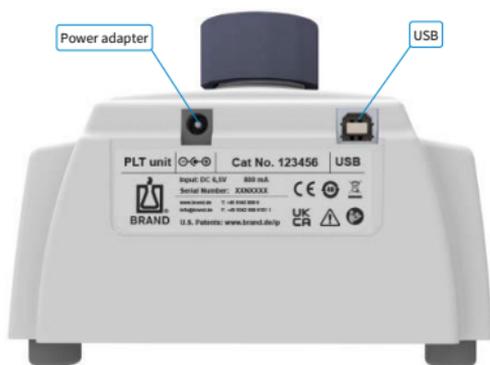

- a. Connect the PLT unit to a mains socket with the mains adapter.
- b. Then connect the PLT unit to a USB 2.0 port on the computer with a connection cable.

4 Operate the program

The PLTconnect program can be started by double-clicking on the BRAND PLTconnect icon on the desktop or via START> All programs> BRAND> PLTconnect .


4.1 Connect to PLT unit

After starting, the program connects automatically with a connected device. If this is successful, “Ready!” is displayed in the lower left status bar.

If the connection to the PLT unit failed, a fault message will be displayed instead.

In order to establish a connection with the device manually, please click on:

File > Connect To PLT unit.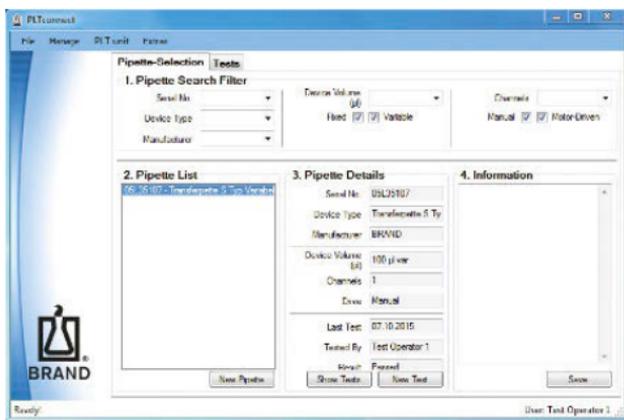


4.2 User settings

The user is shown in the bottom right corner of the user interface.


To create a new user, please click on Manage > Test Operators.

A new tester can be created in the “Manager testers” window or an existing tester can be sought, removed or edited.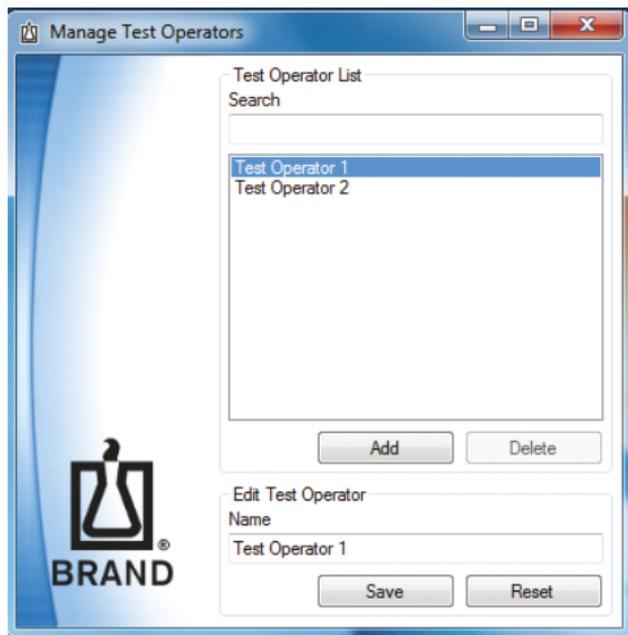

To switch users, please click on File > Switch user.


You can select a user name from the window that opens and confirm this by clicking on Apply.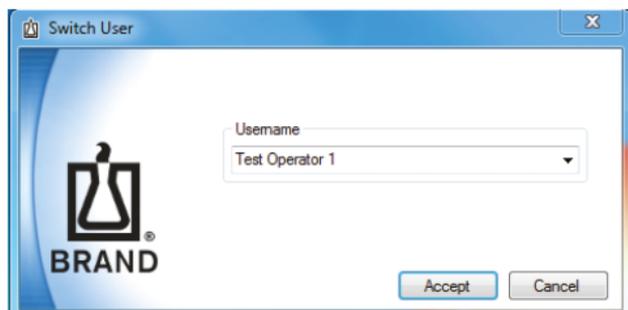

4.3 Creating pipettes


To create a new pipette, please click on the New pipette button on the Pipette selection tab.

Alternatively, you can also select File > Add New Pipette.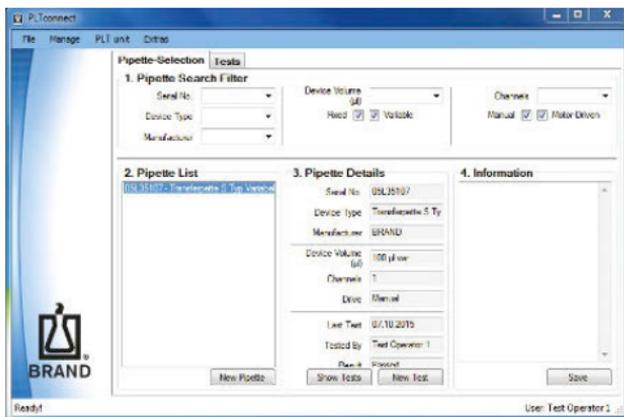

Then follow the instructions from the assistant.

Below: You can filter, insert, remove or change the device types stored in the system under Manage a Device Types.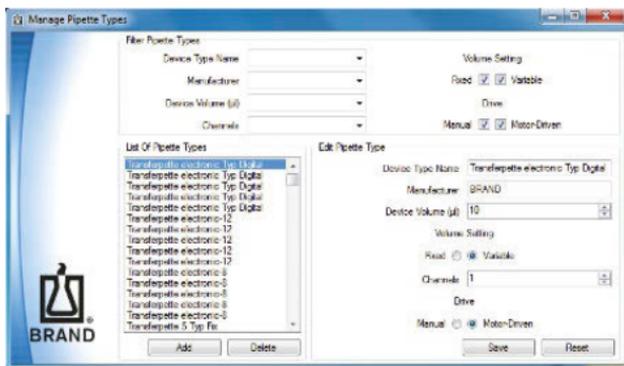


4.4 Carrying out testing

To carry out a test, first click on the New test button on the Pipette selection tab or select File Create new test.

The “Carry out a new test” assistant opens. Please follow the instructions from the assistant.

In the “Test setup – Stipulate test conditions” window, you can determine whether a static or dynamic test is to be carried out and whether it is to be carried out with or without tip installed. The settings will be automatically passed on to the PLT unit.

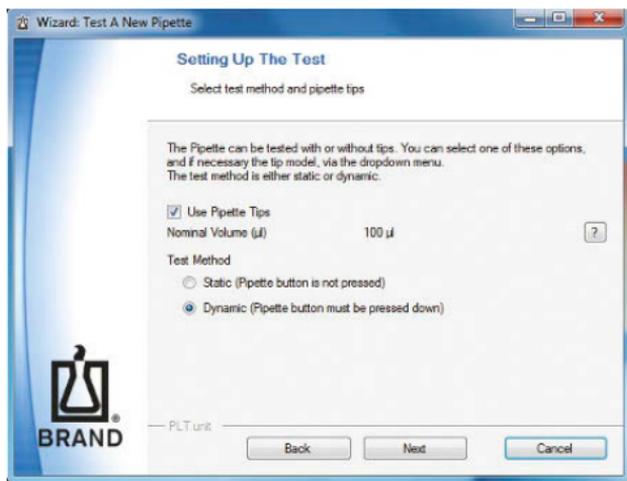
The measurement is started by pressing the left button on the PLT unit.

Then, the result can be transferred to the PC by pressing the right button on the PLT unit.

4 Operate the program

Please tick the “Show Test Certificate” checkbox to display the test certificate. Then you can complete the testing.

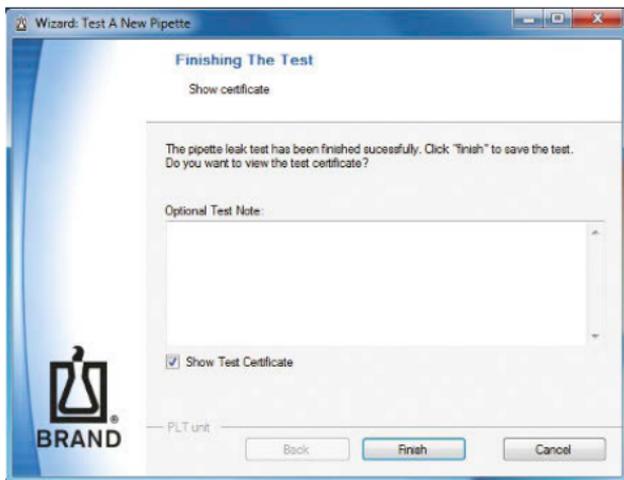BRAND GMBH + CO KG
Postfach/P.O. Box 11 55
D-97877 Wertheim/Main
Germany
<http://www.brand.de>

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
Unterschrift / Signature


4.5 Filter function

You can select pipettes already stored in the system via defined criteria in the “1. Pipette Search Filter” field in the Pipette Selection tab. These are listed in the “2. Pipette List” field.

The details of a pipette selected in field 2 are displayed in the “3. Pipette Details” and in field “4. Information” field. To create or expand a description, enter the desired information into the “4. Information” field and then click on Save.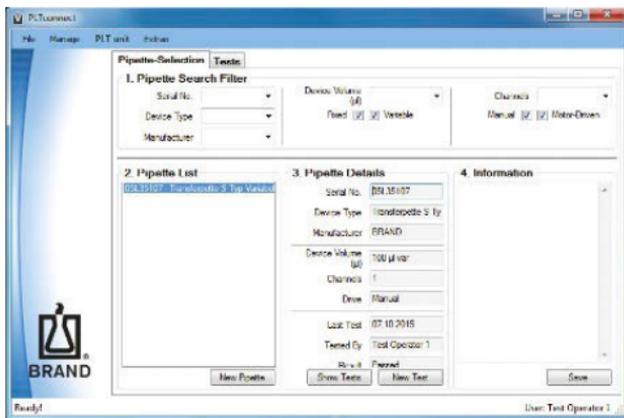

Corresponding field breakdowns and filter functions for displaying tests already carried out can be found in the Tests tab. You can display and print the test certificates here by clicking on the Certificate button.

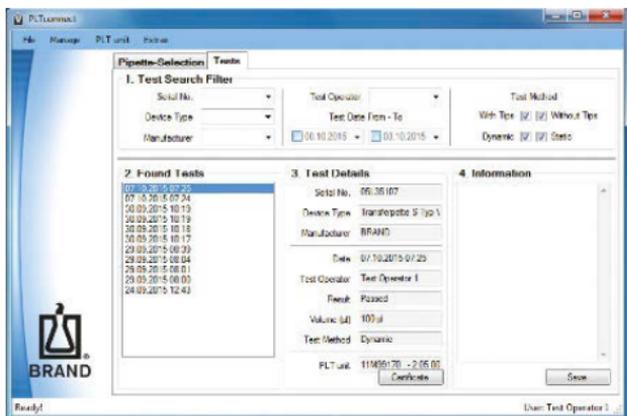4.6 Import/export data

To import existing data sets, please click on:

File > Database Archive > Import From File. and select the corresponding file from the dialog window that opens.

To export datasets, please click on File a Database archive > Export To File.

4.7 Self-test

The self-test is used to detect any leaks in the BRAND PLT unit itself.

There are two different tests available. The quick test takes approx. 2 minutes and should be carried out daily. The full test takes approx. 12 minutes. It includes the additional differentiated testing of the internal tanks and should be carried out weekly.

You can choose between the two tests under PLT unit > Self-test. The appropriate assistant opens.

Follow the instructions from the assistant.

After the test is complete, the test certificate opens.

5 Markings

Symbol or number	Meaning
	With this mark, we confirm that the product complies with the requirements set out in the EC Directives and has been subjected to the specified testing procedures.

6 Disposal

BRAND electronic devices

If you wish to dispose of a BRAND electrical or electronic device, this must be done separately from unsorted municipal waste. This is indicated by the symbol of the crossed-out trash can. BRAND will gladly dispose of your device with this mark.

To do this, please fill out the form for returning old devices at www.brand.de/compliance and send it with the device to the address specified in the form. Please also observe the rules of data protection and, if necessary, delete all data on the devices or USB data carriers before returning them.

| Printed in Germany | 0623-1