HandyStep[®] electronic

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Operating Manual

Please read carefully before use, and follow all operating and safety instructions!

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The device named below fulfills the relevant fundamental requirements of the EC directives and standards listed. This declaration of conformity is issued under the sole responsibility of the manufacturer. In case of unauthorized modifications to the device, the declaration becomes invalid.

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HandyStep® electronic including PD-Tips from BRAND (all sizes) and charging adapter (Identification number 70 50 00 -70 50 04)

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This document declares the accordance with the named harmonized regulations, but does NOT assure specific properties.

目录	页码	Contents	Page	Table des Matières	Page	Contenido	Pág.
安全说明	4	Safety instructions	4	Règles de sécurité	4	Normas de seguridad	4
用途	7	Purpose	7	Utilisation	7	Aplicación	7
使用限制	7	Limitations of use	7	Limites d'emploi	7	Limitaciones de uso	7
禁止操作	8	Operating exclusions	8	Interdictions d'emploi	8	Excepciones de uso	8
操作元件	9	Operating elements	9	Eléments de commande	9	Elementos de manejo	9
初始操作	10	Initial operation	10	Mise en service initiale	10	Puesta en servicio inicial	10
插入电池	10	Insert battery pack	10	Insérer l'unité d'accus.	10	Colocar la unidad de baterías	10
电池充电	11	Charge battery pack	11	Recharger l'unité d'accus.	11	Cargar la unidad de baterías	11
更换电池	14	Replace battery pack	14	Remplacer l'unité d'accus.	14	Reemplazar la unidad de baterías	14
开机/关机	15	Switching on/off	15	Marche et arrêt	15	Encendido y apagado	15
插入吸头	16	Inserting tips	16	Logement des pointes	16	Colocación de las puntas	16
取下吸头	18	Removing tips	18	Enlèvement des pointes	18	Extracción de las puntas	18
HandyStep® electronic		Functions of the HandyStep® electronic	19	Fonctions du HandyStep® electronic	19	Funciones del HandyStep® electronic	19
电子连续分液器 的功能	19	Aspirating liquid	20	Aspiration de milieu	20	Aspiración de muestra	20
吸取液体	20	DISP-mode	22	Mode-DISP	22	Modo-DISP	22
DISP (分液) 模式	22	AUTO-DISP-mode	24	Mode-AUTO-DISP	24	Modo-AUTO-DISP	24
AUTO-DISP (自动分液) 模式	24	PIP-mode	26	Mode-PIP	26	Modo-PIP	26
PIP (移液) 模式	26	Adjusting STEP volume	27	Réglage du volume de pas	27	Ajuste del volumen de paso	27
调整STEP (单次排液) 体积	27	Adjusting aspiration and dispensing speed	28	Réglage de la vitesse d'aspiration et de dosage	28	Ajuste de la velocidad de aspiración y de dosificación	28
调整吸液和分液速度	28	Changing mode	30	Changement de mode	30	Cambio de modo	30
更改模式	30	Use of compatible tips from other manufacturers	31	Utilisation de pointes compatibles d'autres fabricants	31	Utilización de puntas compatibles de otros fabricantes	31
兼容吸头的使用	31	Residual stroke	33	Surplus de course	33	Sobreembolada	33
剩余行程	33	Non-contact dispensing	33	Dosage à jet libre	33	Dosificación de chorro libre	33
非接触式分液	33	Default settings	33	Ajustages d'usine	33	Ajuste de fábrica	33
默认设置	33	Troubleshooting	34	Dérangement - que faire?	34	¿Qué hacer en caso de avería?	36
故障诊断	34	Accuracy table	42	Table de précision	42	Tabla de precisión	42
技术数据	42	Technical data	43	Données techniques	43	Datos técnicos	43
订购信息	43	Ordering data	44	Données de commande	44	Informaciones de pedido	44
准确度表	44	Maintenance and cleaning	46	Entretien et nettoyage	46	Limpieza y mantenimiento	46
保养和清洁	46	UV sterilization	46	Désinfection aux rayons UV	46	Esterilización con UV	46
UV 紫外灭菌	46	Return for repair	46	Retour pour réparation	46	Envíos para reparación	46
返修	46	Contat addresses	48	Adresses de contact	48	Direcciones de contacto	48
联系地址	48	Calibration Service	49	Service de calibrage	49	Servicio de calibración	49
校准服务	49	Warranty · CE note	50	Garantie · Remarque CE	50	Garantía · Nota CE	50
丢弃 · CE	50	Disposal	51	Elimination	51	Eliminación	51
担保信息	51						

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Safety instructions

This instrument may sometimes be used with hazardous materials, operations, and equipment. To address all of the safety problems associated with its use in such applications is beyond the scope of this manual. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Every user must read and understand this operating manual prior to using the instrument and observe these instructions during use.
2. Observe general instructions for hazard prevention and safety instructions; e.g. wear protective clothing, goggles and gloves. When working with infectious or other hazardous samples, all appropriate regulations and precautions must be followed.
3. Observe all specifications provided by reagent manufacturers.
4. Never use the instrument in an atmosphere with danger of explosion. Highly flammable liquids must not be pipetted.

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Safety considerations

- continued -

13. The original battery must not be replaced with non rechargeable batteries or rechargeable batteries of other manufacturers.
14. To charge the NiMH battery pack use only the original AC adapter in combination with the original charging dock.
15. The AC adapter has to be protected against moisture and must be used only for this instrument.
16. Dispose of battery pack only when completely discharged, and according to applicable regulations.

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Note:

Attention: Risk of injury!
Do not touch the plunger seat inside the instrument!

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Warning!

Improper use of the instrument or the batteries (short circuit, mechanical damage, overheating, incorrect AC adapter etc.) may, in extreme cases, lead to explosion of the batteries.

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Purpose

The HandyStep® electronic is a microprocessor controlled battery- operated repetitive pipette. The instrument automatically recognizes new PD-Tips from BRAND, which are encoded for tip volume. This allows fastest possible volume selection. If you want to use compatible tips from other manufacturers, you must set the volume manually.

From now on, the term "tips" refers to PD-Tips with size encoding and compatible tips.

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Limitations of use

The instrument is designed for quick and simple repetitive dispensing of samples, observing the following limitations:

- use between +15 °C and +40 °C (59 °F and 104 °F) (instrument and reagents - other temperatures on request)
- vapor pressure up to 500 mbar
- viscosity：
 - 20 mPa s with 50 ml PD-Tips
 - 260 mPa s with 5 ml PD-Tips
 - 977 mPa s with 1.25 ml PD-Tips

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Operating exclusions

During proper operation, the dispensed sample only comes in contact with the tip and not with the HandyStep® electronic.

The user has to ensure the compatibility of the instrument with the intended application.

Never dispense samples which attack polypropylene, polyethylene or polycarbonate (housing). Avoid aggressive fumes (danger of corrosion)!

The device may not be used for oxidizing acids, since they can attack metallic parts and the electronics.

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Note:

The instrument and the tips are not autoclavable (121 °C / 250 °F). See sterile tips on page 45. Pipette tips are disposable. For disinfection you may use commercially available disinfecting solutions as long as they do not attack PC/PBT or damage the display.

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Initial operation

Insert battery pack

1. Open the battery compartment cover and insert battery pack, ensuring that the electrical contacts on the battery pack connect with those of the HandyStep® electronic.
2. Replace the battery compartment cover.

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Warning!

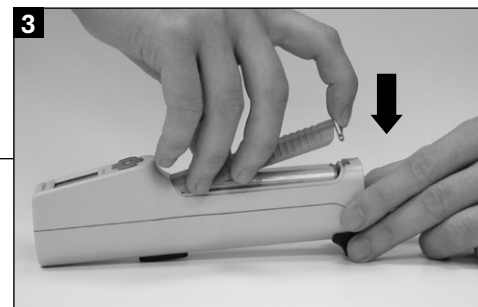
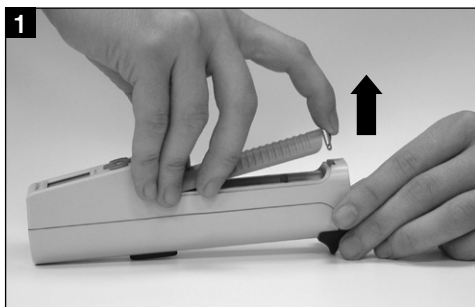
The instrument must not be used or charged in potentially explosive environments.

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Charge battery pack

The battery pack can be charged either inside the HandyStep® electronic or in the charging dock.

Before first use charge the battery pack inside the HandyStep® electronic in the original charging dock for at least 3.5 hours. After 2 - 3 further charging cycles inside the instrument, the full capacity is reached.

Charge indicator is most accurate when battery pack is charged 3.5 hours in the instrument.

Removal of battery pack resets indicator. Upon reinsertion battery-charge frame flashes to indicate that charge status is approximate.

Note:

Stored data and settings are retained when the battery pack is completely discharged or during replacement.

Remove the battery pack when not in regular or daily use.

- A single tone indicates that the battery pack has to be charged or replaced. The battery symbol will flash on the display permanently.

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Charging inside the HandyStep® electronic

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Attention:

Before charging, make sure that the AC adapter is appropriate for your line voltage.

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1. Plug the AC adapter connector into the side of the charging dock.
2. Place the instrument with the battery pack into the charging dock. A tone indicates proper positioning of the instrument.
3. A green light on the charging dock indicates that the instrument is charging.

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The charging time for a nearly empty battery pack mounted in the instrument is approximately 3.5 hours.

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Attention:

Upon each battery pack insertion, charge-indicator frame flashes until battery pack is charged for at least 3.5 hours inside instrument.

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Charging the battery pack in the charging dock

The charging dock allows you to charge a separate battery pack while the HandyStep® electronic is in use.

1. Plug the AC adapter connector into the side of the charging dock.
2. Insert the battery pack into the central opening of the charging dock, making sure the battery pack contacts connect with the charging dock contacts.
3. A green light on the charging dock indicates that the battery pack is charging.

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The charging time for a nearly empty battery pack in the charging dock is approximately 2.5 hours.

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Recharger l'unité d'accus.

– suite –

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绿灯状态	关	开	闪烁
Green light status	Off	On	Flashing
过程	充电基座闲置	正在充电	电池充满; 维持充电
Process	Charging dock empty	Charging in process	Battery pack fully charged; maintenance charge

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Replace battery pack

A fully charged battery pack provides enough power for approximately 2.5 hours of continuous use with a 12.5 ml PD-Tip filled with a sample with similar viscosity and density to water.

- A tone indicates the need for recharging.

- The battery symbol will flash on the display.

This indicates the need to recharge the HandyStep® electronic or to insert a spare battery pack.

Before removing the battery, make sure the tip is completely empty.

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Switching on/off

- The HandyStep® electronic can be switched on by inserting and locking a tip, or by pressing the ENTER key.
- When the locking lever is in the open position, the display shows "CLO" (Close).
- Display: **rEF** (reference run): Press the STEP-key to start reference run.
rEF is displayed after every replacement of the battery pack.

- Auto Power Off

- The HandyStep® electronic is factory-set to switch off after 3 minutes of inactivity.
- To switch off manually, press the ENTER key briefly

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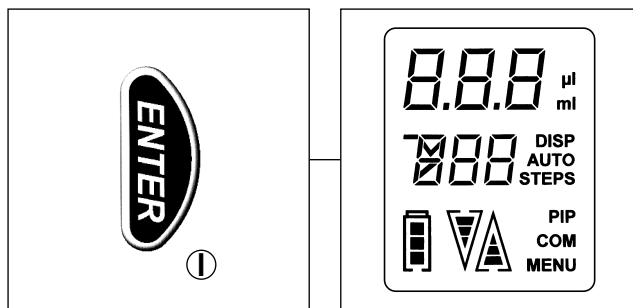
Inserting tips

Attention:

Risk of injury! Do not touch the plunger seat inside the instrument!

1. Push the locking lever to its upper stop.
2. Slide a tip laterally into the Handy-Step® electronic.
3. Push the locking lever down until you hear a distinct clicking (2 steps).
4. If using encoded PD-Tips from BRAND, the PD-Tip capacity and the word **TIP** appear on the display for 3 sec.
5. Then the partial volume to be dispensed in each step is displayed. The default setting is 10 % of the maximum tip capacity.

If using compatible tips (not PD-Tips) you must enter the capacity of the tip yourself. (see page 31)

Note:

When using 25 ml and 50 ml tips, the supplied adapter must be used. It can be connected to the tips using a twist lock and can be removed after use.

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- If the locking lever is not fully engaged, horizontal bars are shown on the display.

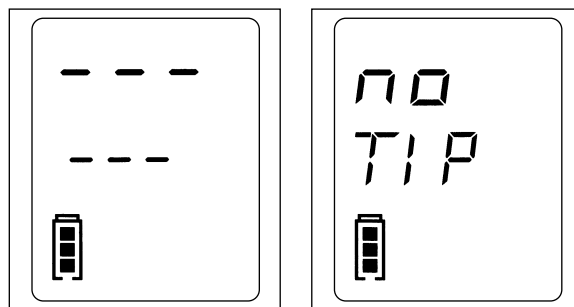
- If the locking lever was fully closed but no tip was inserted, **no TIP** appears on the display.

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Removing tips

1. Hold the HandyStep® electronic over a suitable container.
2. Completely discharge the tip by continuously pressing the STEP key (releasing the residual stroke lock), or by using the "E" mode (see page 23).
3. Push the locking lever to its upper stop.
4. Remove the tip.

Attention:

Don't remove PD-tip before piston of the tip is in the lowest position (tip completely discharged!).

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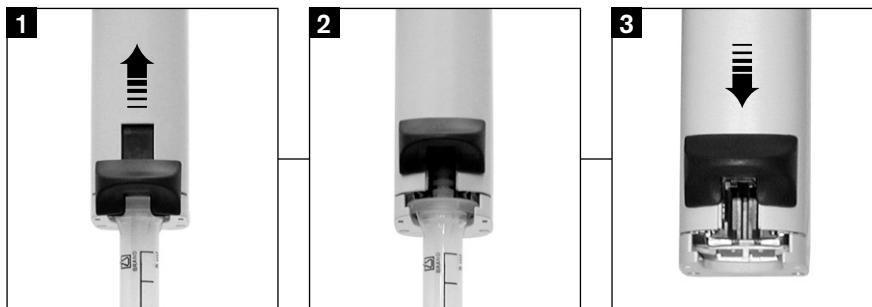
Note:

After dispensing highly viscous or crystallizing sample, the tip may wear and require replacement

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Functions of the HandyStep® electronic

The instrument performs the following functions:

1. Dispensing **DISP** (see page 22):
Default mode after switching on. A user-defined volume can be dispensed repeatedly, depending on the capacity of the tip and the fill-ing volume.
2. Automatic Dispensing **AUTO-DISP** (see page 24):
Identical to DISP-mode, but with automatic dispensing by continuously pressing the STEP key. Releasing the STEP key interrupts the process.
In the AUTO-DISP-mode the instrument has a "Learn" function: After pipetting at least three steps, the HandyStep® electronic calculates the average dispensing interval, and will repeatedly dispense at that interval as long as the STEP key is pressed.
3. Pipetting **PIP** (see page 26):
In PIP-mode, the HandyStep® electronic functions as a positive displacement non-repeating pipette, particularly useful for volatile, foamy or high viscous materials. A preselected volume is aspirated and dispensed (Positive displacement principle).

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Note:

Default setting for the HandyStep® electronic is DISP mode (see page 22).

We recommend that a new PD tip has to be primed before use. For this, fill the tip at a minimum with the liquid and then empty it again (see page 21). Small air bubbles in the area of the piston after priming do not affect the results.

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1. The arrow on the display points upward △ ("Aspirate").
2. Push the STEP key once. The tip will be filled completely.
3. During filling, the number of remaining steps is continuously displayed.
4. After the filling procedure has been completed, the arrow on the display shows ▽ ("Dispense").

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To interrupt the filling procedure and to bleed the PD tip:

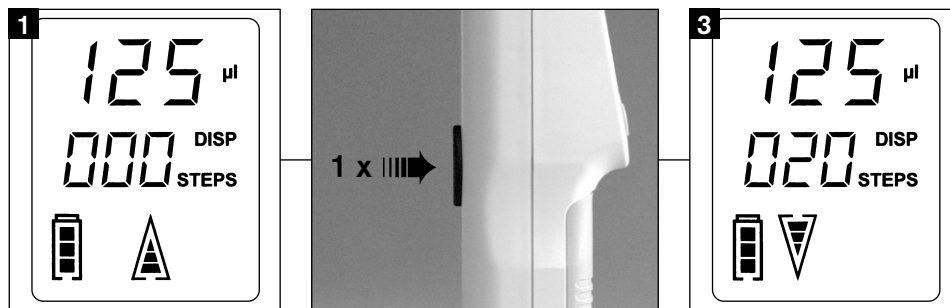
To interrupt the filling process, press the STEP key. To continue filling, press the Δ key first and hold down the STEP key in succession. The step count on the display will increase. Activate the ∇ button and press and hold down the STEP button to completely discharge the liquid again.

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Attention:

After aspirating liquid, the Handy-Step® electronic will automatically dispense a small amount of liquid in order to compensate for gear backlash.

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1. The arrow on the display points downward ▽ ("Dispense").
2. Press the STEP key to dispense the selected step volume (see "Adjusting STEP volume" page 27).
3. The display shows the remaining number of steps.
4. After dispensing the last full step, a small amount of sample will be left in the tip. To discharge this volume, hold down the STEP key.
5. Remove tip (see page 18).

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Note:

The tip must be completely empty before you remove it.

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Attention:

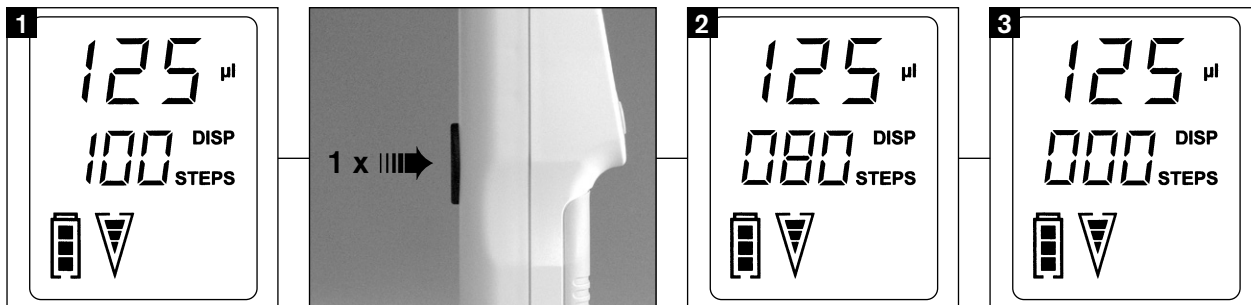
Dispensing can be interrupted at any time by pressing the △ key.

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Note:

To empty the tip at any time, press the ▽ key and then hold down the STEP key. The letter E appears on the display and the arrow pointing downwards flashes. This process can be interrupted at any time by pressing the △ key.

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A partially filled tip can be completely filled by pressing and holding, in sequence, the △ key and the STEP key.

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AUTO-DISP-mode

1. To activate the AUTO-DISP-mode, briefly press the MENU key twice.

Note:

The AUTO-DISP-mode must be actively selected.

2. Select the AUTO-DISP-mode with the ▽△ keys (ON/OFF).
3. Confirm by pressing ENTER.
4. "Learn" feature: After pipetting at least three steps, the HandyStep® electronic automatically determines the average time interval between steps (page 19).
5. Hold down the STEP key to dispense the selected volume (page 27) at the calculated intervals.

Note:

Intervals can range from 0.1 s to 5 s.

6. To interrupt dispensing, release the STEP key. The calculated step interval remains in memory.
7. In AUTO-DISP-mode, the "Learn" process can be restarted at any time.

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PIP-mode

(Reverse pipetting,
positive displacement principle)

1. Select the PIP-mode by holding down the MENU key (> 3s), and press it briefly (page 30).
Make sure that plunger is at lowest position!
2. Select a volume (page 27).
3. To aspirate the selected volume, press the STEP key. A volume slightly larger than the selected volume will be drawn into the tip.
4. To dispense the volume, press the STEP key again. Some residual volume will remain in the tip.
5. To aspirate a new sample, press the STEP key again.
6. When you have completed pipetting, discharge the residual stroke by holding down the STEP key.

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Attention!

To interrupt the pipetting process at any time, press the Δ key.

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Note:

To change from PIP-mode to AUTO-DISP-mode, you have to switch to DISP-mode first.

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Adjusting STEP volume

Note:

After the tip size has been recognized, the dispensing volume is shown on the display.

The default step volume is 10 % of the nominal capacity.

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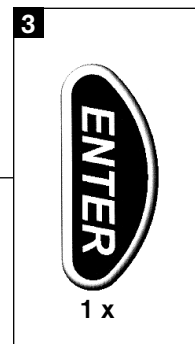
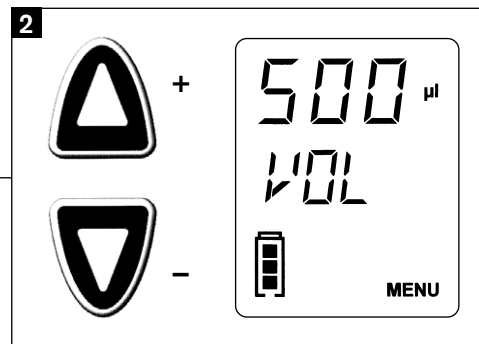
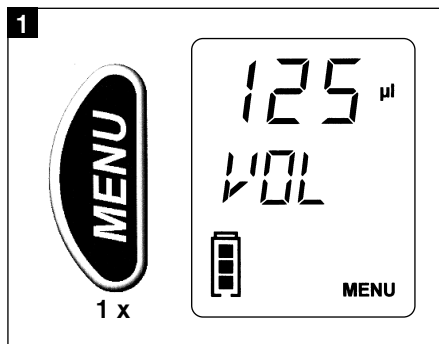
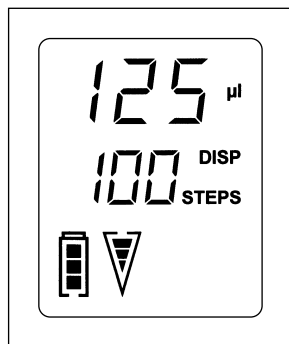
1. Press the MENU key once. The volume (VOL) is shown on the display.
2. Use the ▽△ keys to select the appropriate volume. Holding down the ▽△ keys will accelerate the rate of change.
3. Confirm the volume selection by pressing ENTER.

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Adjusting aspiration and dispensing speed

1. Hold down the MENU key (> 3s) to adjust the aspiration speed. The arrow for selection of aspiration speed lights up on the display. Choose between three different aspiration speeds.

Note:

Lowest speed:

Display shows outline of arrow with one bar.

Highest speed:

Display shows outline of arrow with three bars.

2. Use the $\nabla\Delta$ keys to select aspiration speed.
3. Confirm aspiration speed setting by pressing ENTER.
4. The down arrow appears on the display.
5. Use the $\nabla\Delta$ keys to select dispensing speed.
6. Confirm settings by pressing ENTER. You are returned to the main menu.

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Use of compatible tips

Compatible dispenser tips which are not manufactured by BRAND are not encoded for size. These tips can still be used with the HandyStep® electronic (see "Insert tips" page 16). After insertion of one of these tips, the display will automatically go to tip selection mode.

By default, the 5 ml tip is displayed flashing regardless of the inserted tip. 12 possible tip sizes are available.

1. Use the ▽△ keys to select appropriate tip size.
2. Confirm setting by pressing ENTER.

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– continued –

Attention!

For convenience, when inserting the next unencoded tip, the previously selected tip size will come up first flashing on the menu. Confirm by pressing ENTER.

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At the time of printing of this operating manual, the following dispenser tips were compatible with the HandyStep® electronic:

- PD-Tips
(BRAND GMBH + CO KG, Germany)
- Encode™ tips
(Rainin Instrument LLC, USA)
- Repet-Tips
(Gilson S.A.S., France)
- Combitips®, Combitips® plus, Combitips advanced®
(Eppendorf AG, Germany)
- ritips®
(ritter GmbH, Germany)
- Distritips®
(Gilson S.A.S., France)
- Fisherbrand® Dispenser-Tips
(Fisher Scientific, USA)
- VWRbrand® Combi-Syringes
(VWR Scientific Holding Corp., USA)

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Residual stroke

A lock prevents the residual stroke from being dispensed as a regular dispensed or pipetted volume. The symbol Δ indicates this remaining volume. It can be discharged by holding the STEP key.

Non-contact dispensing

When tips with volumes of 5 ml or larger are selected, the default dispensing speed adjusts to allow accurate dispensing without contacting the walls of vessel (free-jet dispensing). This can be overridden by selecting a different speed (page 28).

Default settings

To revert the HandyStep® electronic to its factory default settings, simultaneously press the MENU and ENTER keys for > 3 seconds, with the plunger in the lowest position.

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Troubleshooting

Problem	Error message displayed	Possible cause	Corrective action
Tip dripping	No message	Leaking tip	Replace tip
Faint display; reduced speed	Battery symbol flashing	Battery low	Recharge battery pack or replace with charged battery pack.
Instrument not responding	ERR 001	- Locking lever not engaged during aspirating or dispensing - Locking lever open at lower third of the tip	1. Press ENTER to confirm error. Follow MENU-instructions to avoid instrument damage! Display: CLO (Close) 2. Engage the locking lever Display: rEF (Reference run) 3. Press the STEP key Attention: Liquid remaining in tip will be discharged during the reference run (rEF). Discharge sample into a suitable vessel (see pages 5, point 10).
Instrument not responding	ERR 002	Motor blocked - Battery low - Viscosity is too high (see page 7) - Tip orifice obstructed	1. Press ENTER to confirm error. Follow MENU-instructions to avoid instrument damage! - Recharge or replace battery pack - Reduce aspirating/dispensing speed Display: rEF (Reference run) 2. Press the STEP key Attention: Liquid remaining in tip will be discharged during the reference run (rEF). Discharge sample into a suitable vessel (see pages 5, point 10).
Instrument not responding	ERR TIP	Locking lever opened too early within the lower third of the tip	1. Press ENTER to confirm error. Follow MENU-instructions to avoid instrument damage! 2. Display: OPE (Open) Push the locking lever to its upper stop.

Troubleshooting - continued -

Problem	Error message displayed	Possible cause	Corrective action
Instrument not responding - continued -			3. Display alternating between out TIP and CLO . Remove tip and close locking lever (without tip). 4. HandyStep® electronic automatically carries out a reference run. 5. Display: no TIP Insert new tip (see page 16).
Instrument not responding	ERR 003	Electronic fault or tip connector not at lowest position when inserting new tip	1. Press ENTER to confirm error. 2. Follow MENU-instructions: - Display rEF (reference run) - Press STEP-Key Attention: Liquid remaining in tip will be discharged during reference run (rEF). Discharge sample into suitable vessel (see page 5, point 10). If error message is still shown on the display remove battery pack and insert it again (reset). If this does not solve the problem send instrument back for repair.
No tone, no light when placing the HandyStep® electronic into the charging dock	ACC	- Instrument contains no battery - Battery pack inserted improperly - Battery pack faulty	- Insert battery pack. - Make sure battery contacts are aligned with contacts in the instrument. - Replace battery pack.
Charge-indicator is flashing	No message	Removal or insertion of a new battery pack	- Charge battery pack for at least 3.5 hours inside the HandyStep® electronic.
Dispensing process interrupted. Attention! Volume not completely discharged	ERR VOL	Dispensing process interrupted by - Pressing the Δ key during dispensing - electronic faults	1. Press ENTER to confirm error. 2. Restart dispensing by pressing the STEP key.

Note:

If other error messages are displayed, remove battery pack and insert it again.
 If this does not solve the problem, send instrument in for repair.

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Accuracy table

Measurements were carried out using original BRAND PD-Tips and distilled water.

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The nominal volume is the maximum volume printed on the PD-Tip.

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The predetermined ISO 8655 tolerances were not exceeded.

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Note:

The device is marked in accordance with the German Measurement and Calibration Act as well as the Measurement and Calibration Regulation:

DE-M 20

Character string: DE-M (DE for Germany), framed by a rectangle, as well as the last two digits of the year in which the marking was affixed (here: 2020).

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Technical data

Battery pack:

Rechargeable Nickel Metal Hydride battery pack:
NiMH, 4 x AAA cells,
4.8 V/700 mAh

Charging times for empty battery pack:

in battery compartment of charging dock: approx. 2.5 h

inside the HandyStep® electronic: approx. 3.5 h

When battery capacity starts to diminish, recharge battery pack in the battery compartment of the charging dock.

Battery pack capacity:

A full charge gives approx. 2.5 hours of continuous work with a 12.5 ml PD-Tip.

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Maintenance and cleaning

The HandyStep® electronic is calibrated at the factory and is maintenance-free. If required, clean the instrument with a cloth dampened with water or a mild soap solution. **The HandyStep® electronic must not be dismantled.**

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UV sterilization

The unit can withstand the usual output of a UV sterilization lamp. The effects of the UV may cause some color change.

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Return for Repair

Attention:

Transporting of hazardous materials without a permit is a violation of federal law.

- Clean and decontaminate the instrument carefully.
- It is essential always to include an exact description of the type of malfunction and the media used. If information regarding media used is missing, the instrument cannot be repaired.
- Shipment is at the risk and the cost of the sender.

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– continued –

Outside the U.S. and Canada:

- Complete the “Declaration on Absence of Health Hazards” and send the instrument to the manufacturer or supplier. Ask your supplier or manufacturer for the form. The form can also be downloaded from www.brand.de.

In the U.S. and Canada:

- Contact BrandTech Scientific, Inc. and obtain authorization for the return **before** sending your instrument for service.
- Return only cleaned and decontaminated instruments, with the Return Authorization Number prominently displayed on the outside of the package to the address provided with the Return Authorization Number.

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India:

BRAND Scientific Equipment
Pvt. Ltd.
303, 3rd Floor, 'C' Wing, Delphi
Hiranandani Business Park, Powai
Mumbai - 400 076 (India)
Tel.: +91 22 42957790
Fax: +91 22 42957791
E-Mail: info@brand.co.in
www.brand.co.in

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Calibration Service

ISO 9001 and GLP-guidelines require regular examinations of your volumetric instruments. We recommend to check the volume every 3-12 months. The interval depends on the specific requirements on the instrument. For instruments frequently used or in use with aggressive media, the interval should be shorter.

The detailed testing instruction can be downloaded on www.brand.de or www.brandtech.com.

BRAND also offers you the possibility, to have your instruments calibrated by the BRAND calibrating service or the BRAND-owned DAkKS calibration service.

Just send in the instruments to be calibrated accompanied by an indication of which kind of calibration you wish. Your instruments will be returned within a few days together with a test report (BRAND calibrating service) or with a DAkKS calibration certificate. For further information please contact your dealer or BRAND.

Complete ordering information is available for download at www.brand.de (see technical documentation).

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Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operation or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual.

We are not liable for damage resulting from any actions not described in the operating manual or if non-original spare parts or components have been used.

U.S. and Canada:

Information for warranty please see www.brandtech.com.

Subject to technical modification without notice. Errors excepted.

This sign certifies that the product meets the requirements of the EEC directive and has been tested according to the specified test methods.

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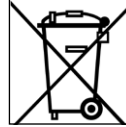
Disposal

The adjoining symbol means that storage batteries and electronic devices must be disposed of separately from household trash (mixed municipal waste) at the end of their service life.

- According to the Directive 2002/96/EC of the European Parliament and of the Council on Waste Electrical and Electronic Equipment (WEEE) of 27 January 2003, electronic equipment requires disposal according to the relevant national disposal regulations.
- Batteries contain substances that can have harmful effects on the environment and human health. Therefore according to the Directive 2006/66/EC of the European Parliament and the Council on Waste Batteries of 6 September 2006 batteries require disposal according to the relevant national disposal regulations. Dispose of batteries only when completely discharged.

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Warning!

Do not short-circuit the battery to discharge it!

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

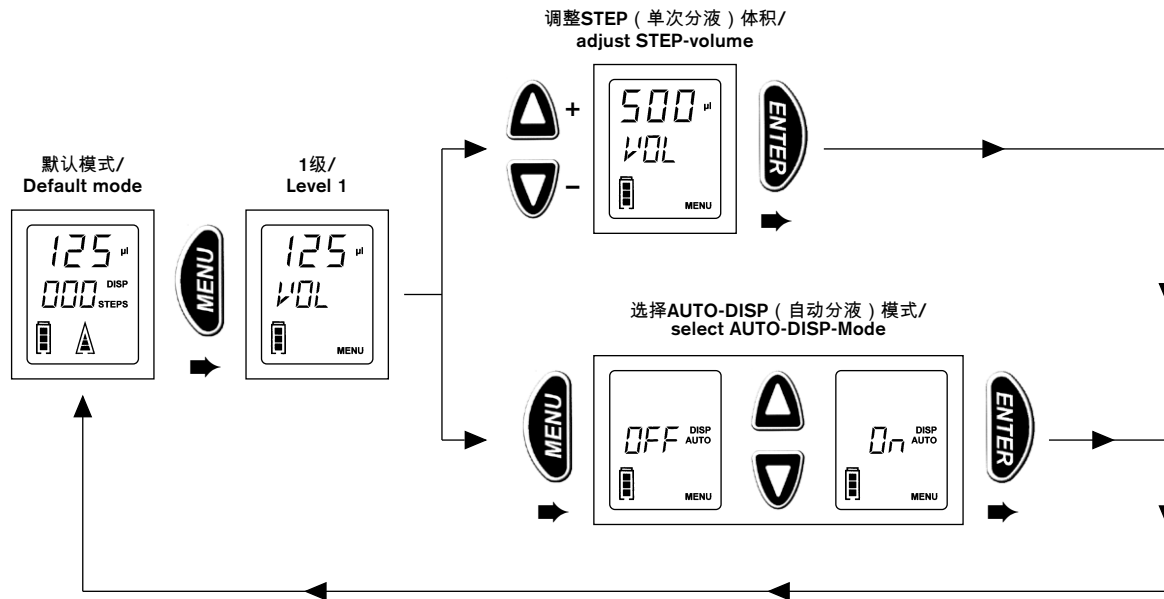
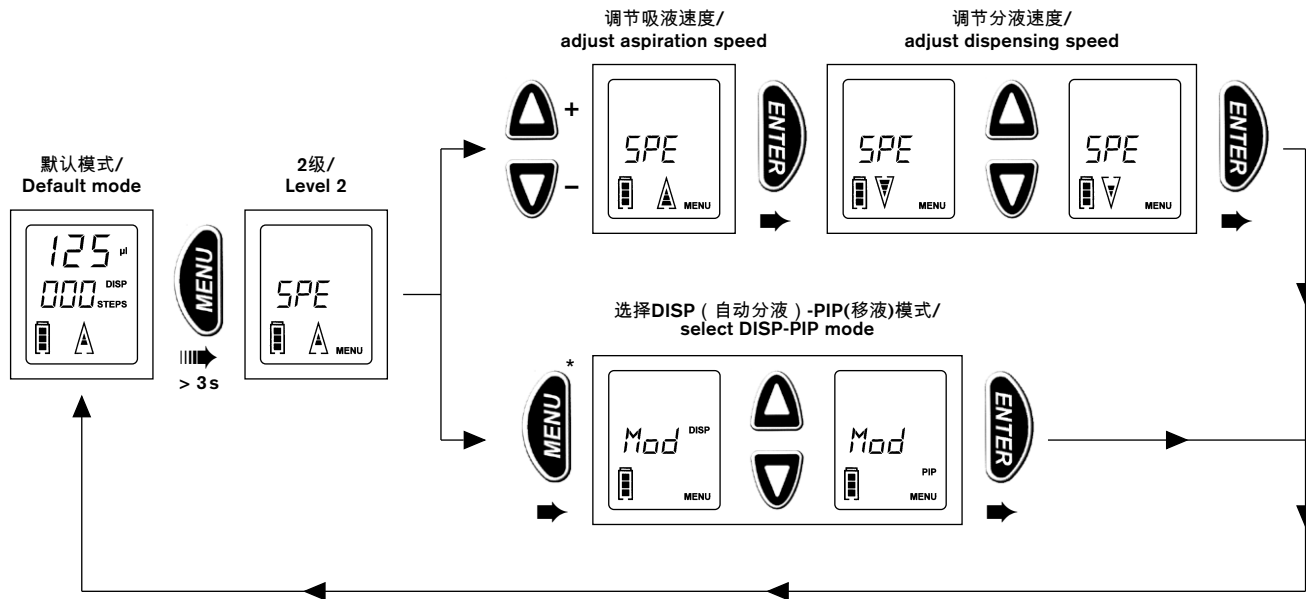
* To select this mode the PD-Tip must be completely empty!

- ➡ Appuyer brièvement sur le bouton MENU / Accionar el botón MENU
- ⏏ Maintenir enfoncé le bouton MENU (> 3s) / Accionar el botón MENU continuellement (> 3s)

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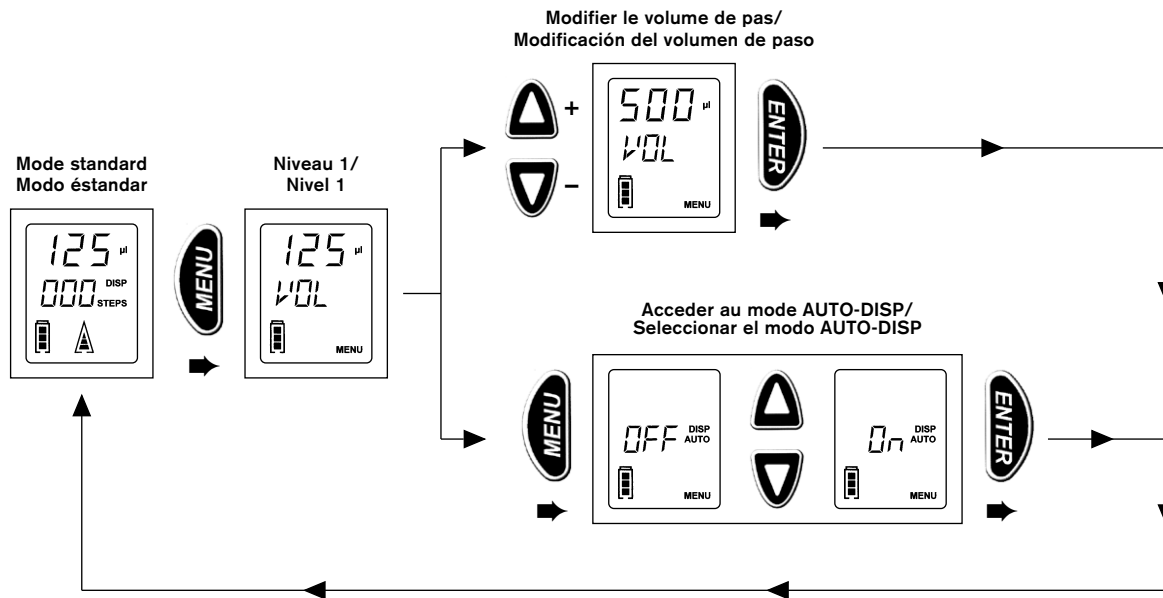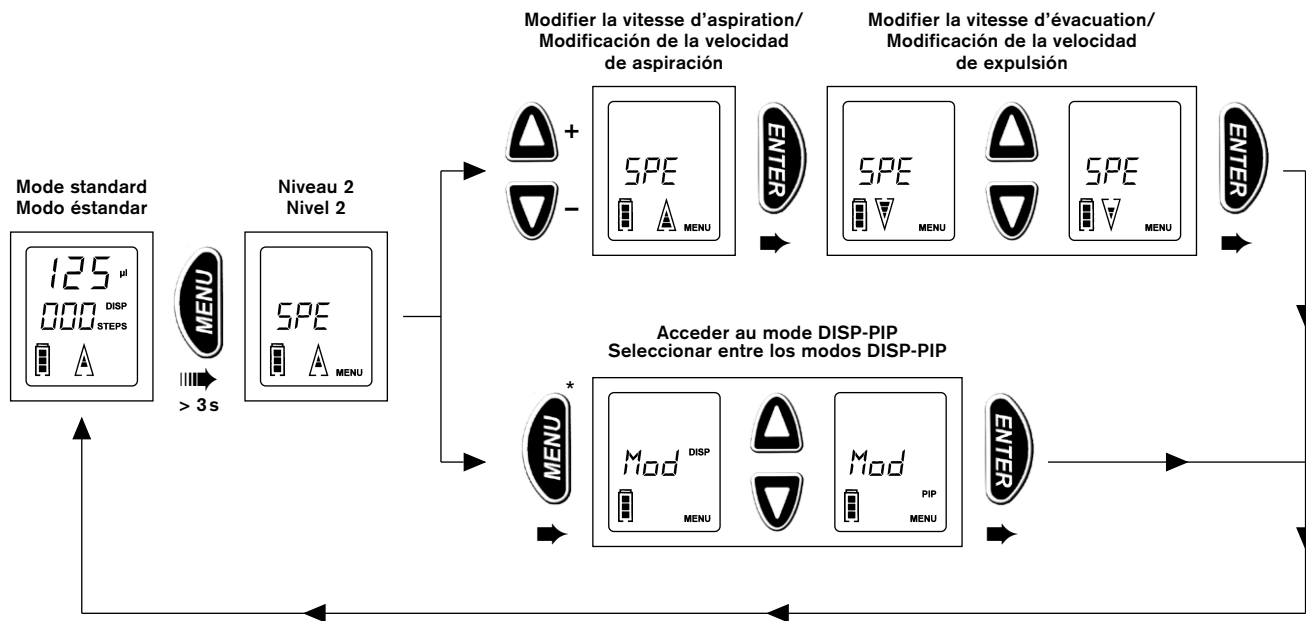BRAND GMBH + CO KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in BRAND products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a "Product Conformity Assessment" (PCA) procedure was performed. As defined in GB/T 26572 the "Maximum Concentration Value" limits (MCV) apply to these restricted substances:

- | | |
|---|-------|
| ● Lead (Pb): | 0.1% |
| ● Mercury (Hg): | 0.1% |
| ● Cadmium (Cd): | 0.01% |
| ● Hexavalent chromium (Cr(+VI)): | 0.1% |
| ● Polybrominated biphenyls (PBB): | 0.1% |
| ● Polybrominated diphenyl ether (PBDE): | 0.1% |

Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The environmental Friendly Use Period for BRAND instruments is 40 years.

此表格是按照SJ/T 11364-2014中规定所制定的。
This table is created according to SJ/T 11364-2014.

MATERIAL CONTENT DECLARATION FOR BRAND PRODUCTS							
有毒有害物质或元素 Hazardous substances							
部件名称 Part name	铅 Pb	汞 Hg	镉 Cd	六价铬 Cr(+VI)	多溴联苯 PBB	多溴二苯醚 PBDE	环保期限标识 EFUP
包装 / Packaging	O	O	O	O	O	O	
塑料外壳 / 组件 Plastic housing / parts	O	O	O	O	O	O	
电池 / Battery	O	O	O	O	O	O	
玻璃 / Glass	O	O	O	O	O	O	
电子电气组件 Electrical and electronic parts	X	X	X	O	O	O	
金属外壳 / 组件 Metal housing / parts	X	O	O	O	O	O	
电机 / Motor	X	O	O	O	O	O	
配件 / Accessories	X	O	O	O	O	O	

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Note: Table applies to all products. Some of the components or parts listed above may not be part of the enclosed product.

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
- O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
- X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

除上表所示信息外，还需声明的是，这些部件并非是有意图用铅 (Pb) 汞 (Hg)，镉 (Cd)，六价铬 (Cr(+VI))，多溴联苯 (PBB) 或多溴二苯醚 (PBDE) 来制造的。
Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (Cr(+VI)), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by BRAND may enter into further devices or can be used together with other appliances. With these third party products and appliances in particular, please note the EFUP labeled on these products. BRAND will not take responsibility for the EFUP of those products and appliances.

Place, date: Wertheim, 25/02/2019

Hans-Walter Kern (Managing Director Logistics and Production)	i. A. Josef Frohl (Quality Management)
--	---

BRAND GMBH + CO KG · Otto-Schoft-Str. 25 · 97877 Wertheim · Germany
Tel.: +49 9342 808-0 · Fax: +49 9342 808-98000 · E-Mail: info@brand.de
Internet: www.brand.de

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