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Before initial operation:

Observe the Operating Manual, particularly safety instructions, exclusions and limitations of the Dispensette® S model used.

Attention!

The flexible discharge tubing must not be used with hydrofluoric acid (HF).

Operating Exclusions:




Flexible discharge tubing for Dispensette® S must not be used with:

- liquids attacking borosilicate glass (tube holder) or fluoroplastics like PFA and PTFE (e.g., dissolved sodium azide*)
- Hydrofluoric acid
- trifluoroacetic acid
- explosive liquids (e.g., carbon disulfide)
- suspensions (e.g., of charcoal) as solid particles may clog or damage the instrument

* Dissolved sodium azide permitted up to a concentration of max. 0.1%.

Attention!

There should be no visible damage to the discharge tube (e.g. kinks or the like). Each time you are going to use the tubing, examine it carefully! To dispense aggressive liquids, you should take safety measures in addition to the normal precautions. We recommend the use of a protective shield. For cleaning rinse the tube carefully. Do not dismantle!

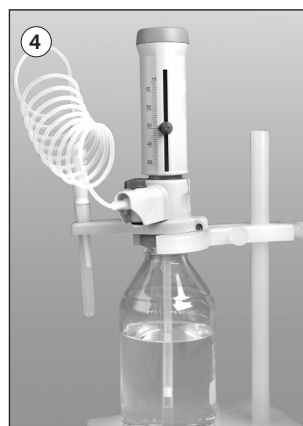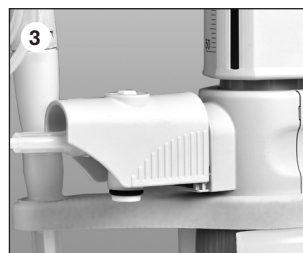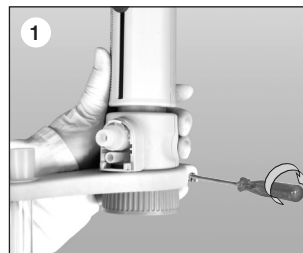Assembly:

1. If the Dispensette® S was already in use, the instrument must be cleaned before mounting the flexible discharge tube.
2. Instruments with a recirculating valve should be set to 'Recirculate', and the valve lever pulled upwards to remove it.
3. Slide the discharge tube housing all the way up. Then pull the discharge tube forward with gentle up and down motions.
4. Push the flexible discharge tube holder from the bottom of the valve block (Fig. 1) and tighten it. For this, Dispensette® S must not be mounted on the bottle. Install the receiver tube.
5. Press the cock plug of the recirculation valve downwards (Fig. 2).
6. Slide the flexible discharge tube housing into the valve block up to the stop (Fig. 3).
7. Slide the discharge tube housing all the way down.
8. Attach the valve lever and press in firmly (Fig. 4). Note the color coding and valve marking of the valve lever (see table below).

Discharge tube	Valve lever
without valve marking	red
ORG	yellow
Pt-Ir	grey – Pt-Ir
TA	grey – Ta

Note:

The coil of the tubing can be stretched to a length of the 800 mm max. The entire coil must lie in regular loops and must not be twisted.

When working with the flexible discharge tube, the specified accuracy and coefficient of variation of the instrument are only obtained for volumes > 2 ml and by gently approaching the upper and lower stops.

Attention!

The bottle must be supported using a bottle stand to help prevent the bottle from accidentally tipping over. To help avoid reagent splashing from the tube, always grip the tube firmly by the handle and replace in the holder after use.

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