

For optimal fluorescence and spectral photometric analysis in the UV/Vis range




Cuvette

4 clear sided

F I R S T C L A S S · B R A N D

Plastic disposable UV-Cuvettes for the UV/Vis range. UV-transparent plastic cuvettes replace fragile glass or quartz cuvettes in many applications. Designed for single use, they eliminate time wasting washing, and the risk of cross-contamination.

- Four optical windows for precise measurements
- Suitable for fluorescence spectroscopy
- Significantly lower costs for use when compared to quartz glass cuvettes
- UV-Cuvettes show minimal autofluorescence
- Will fit most fluorescence photometers
- UV-Cuvettes show very high chemical resistance
- UV-Cuvettes for measurements from 230 nm
- Frosted end for labelling

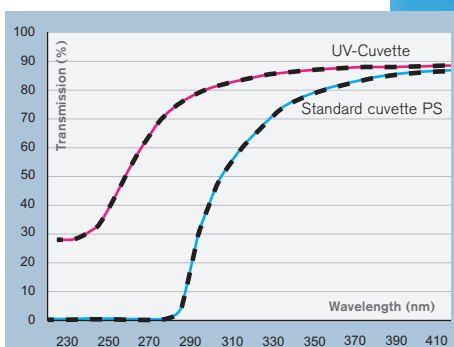

Technical Information

Filling volume:

PS Cuvette: 2.5 - 4.5 ml

UV-Cuvette: 2.5 - 4.5 ml

Dimensions:


Window (W x H): 10 x 35 mm

Light path: 10 mm

Wavelength:

PS Cuvette: 340-900 nm

UV-Cuvette: 230-900 nm

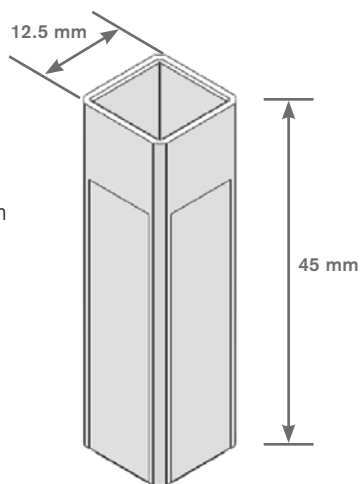

Ordering Data

4 clear sided cuvettes

Description	Pack of	Cat. No.
Cuvette, macro (PS), 4 clear sided	100	7590 30
	500	7590 35
UV-Cuvette, macro, 4 clear sided	100	7591 25
	500	7591 28

Cuvette rack

PP, gray. Holds 16 cuvettes, numbered positions.

L x W x H: 210 x 70 x 38 mm.

Autoclavable (121 °C).

Pack of 1.

Cat. No.	7595 00
----------	---------

Overview Table

Cuvette type	Filling volume min.	Filling volume max.	Dimensions window (W x H)	Range of application	Standard deviation in extinction units
UV-Cuvette, macro, 4 clear sided	2.5 ml	4.5 ml	10 x 35 mm	230 to 900 nm	240 nm $\leq \pm 0.007$ 300 nm $\leq \pm 0.005$
Cuvette, macro (PS) 4 clear sided	2.5 ml	4.5 ml	10 x 35 mm	340 to 900 nm	360 nm $\leq \pm 0.005$


BRAND also manufactures standard macro and semi-micro cuvettes in PS and PMMA. Detailed information on request, or at www.brand.de

Chemical resistance* of plastic cuvettes

Substance	PS	PMMA	UV-Cuvette
Acetic acid, 100%	-	-	+
Acetone	-	-	+
Ammonia	+	+	+
Benzaldehyde	-	-	+
Butanone	-	-	+
Chloroform	-	-	-
Dioxane	-	-	+
DMF	-	-	+
DMSO	-	-	+
Ethyl acetate	-	-	+
Hexane	-	+	-
Hydrochloric acid, 36%	+	-	+
Hydrofluoric acid, 10%	+	+	+
Isopropanol	+	+	+
Nitric acid, 65%	-	-	+
Sodium hydroxide	+	+	+

* Short time resistance, 30 min. Longer term storage of these chemicals should be confirmed by the user. Free samples upon request.

UV-Cuvettes of minimal autofluoreszenz

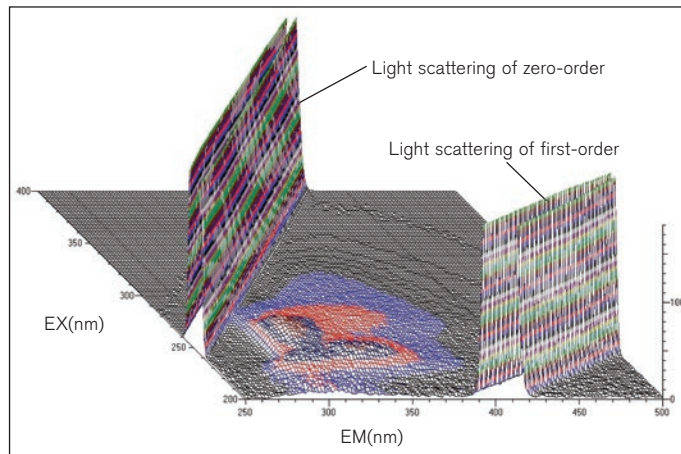

3-D scan from 200 to 400 nm wavelength with Hitachi F-7000 FL-Spectrometer

BRAND®, BRAND. For lab. For life.® and the BRAND word and figurative mark are registered trademarks of

BRAND GMBH + CO KG, Germany.

Our technical literature is intended to inform and advise our customers. However, the validity of general empirical values, and of results obtained under test conditions, for specific applications, depends on many factors beyond our control. Please appreciate, therefore, that no claims can be derived from our advice. The user is responsible for checking the appropriateness of the product for any particular application.

California Residents: For more information concerning California Proposition 65, please refer to www.brand.de/calprop65

Subject to technical modification without notice. Errors excepted.

BRAND GMBH + CO KG · P.O. Box 11 55 · 97861 Wertheim · Germany

Tel: +49 9342 808-0 · Fax: +49 9342 808-98000 · E-Mail: info@brand.de · Internet: www.brand.de

