

Relax – everything is under control!

Pipetting has never been so comfortable

Comfortable

Ergonomic handgrip, weight only 190 g, perfectly balanced design – for fatigue-free pipetting even in prolonged operations.

Sensitive

The *accu-jet® pro* offers motor speed control to set the range of pipetting speeds; use the sensitive pipetting buttons for continuously variable pipetting speed within the selected range.

Powerful

At maximum motor speed, a 50 ml pipette is filled in less than 10 seconds. Now that is fast!

Quiet

Motor and pump operate with quiet discretion. The longer you use it, the more you will appreciate it.



Pipetting keys

Recharging-indicator (LED)

Battery compartment cover

Mode selection
(gravity-delivery/blow-out)

Variable motor speed

Dispensing liquid

Choose between gravity delivery (for pipettes calibrated "to deliver") or blow-out delivery with motor power.

Recharging-indicator

The flashing recharging-indicator LED alerts you that the battery is running low. Don't worry – you can work for another two hours before it's empty – you can use the instrument while it's charging.

Motor speed

Select the desired maximum pipetting speed by turning the thumbwheel. A slower speed setting provides sensitive, controlled operation with lower volume pipettes.

Battery compartment cover

The *accu-jet® pro* is supplied with two battery compartment covers. The second cover enables you to place the instrument inverted on the bench, to avoid contamination of the pipette adapter.