

Dispenser selection chart

| Reagent | seripettor® | seripettor® pro |
|-----------------------------|-------------|--------------------|
| Acetaldehyde | | + |
| Acetic acid, 5% | + | + |
| Acetic acid, 96% | | + |
| Acetic acid (glacial), 100% | | + |
| Acetone | | + |
| Acetonitrile | | + |
| Acetophenone | + | |
| Acetylacetone | + | + |
| Acrylic acid | | + |
| Acrylonitrile | | + |
| Adipic acid | + | + |
| Agar (60 °C) | + | |
| Allyl alcohol | + | + |
| Aluminium chloride | + | + |
| Amino acids | + | + |
| Ammonia, 30% | + | + |
| Ammonium chloride | + | + |
| Ammonium fluoride | + | + |
| Ammonium sulfate | + | + |
| Amyl alcohol (Pentanol) | + | + |
| n-Amyl acetate | | + |
| Aniline | | + |
| Barium chloride | + | + |
| Benzaldehyde | | + |
| Benzyl alcohol | | + |
| Benzylamine | | + |
| Benzylchloride | | + |
| Boric acid, 10% | + | + |
| BSA serum | + | + |
| Butanediol | + | + |
| 1-Butanol | | + |
| Butylamine | | + |
| n-Butyl acetate | | + |
| Calcium carbonate | + | + |
| Calcium chloride | + | + |
| Calcium hydroxide | + | + |
| Calcium hypochlorite | | + |
| Chloroacetaldehyde, 45% | | + |
| Chloroacetic acid | | + |
| Chromic acid, 50% | | + |

| Reagent | seripettor® | seripettor® pro |
|--------------------------------|-------------|--------------------|
| Copper sulfate | + | + |
| Cumene (Isopropyl benzene) | | + |
| Diethylene glycol | + | + |
| Dimethyl sulfoxide (DMSO) | | + |
| Dimethylaniline | | + |
| Ethanol | + | + |
| Ethidium bromide solution | + | + |
| Formaldehyde, 40% | + | + |
| Formamide | + | + |
| Formic acid, 100% | | + |
| Glycerol | + | + |
| Glycol (Ethylene glycol) | + | + |
| Glycolic acid, 50% | + | + |
| Guanidine hydrochloride | + | + |
| HEPES buffer | + | + |
| Hexanoic acid | + | + |
| Hexanol | | + |
| Hydriodic acid | + | + |
| Hydrobromic acid | | + |
| Hydrochloric acid, 37% | | + |
| Hydrogen peroxide, 35% | + | |
| Isoamyl alcohol | | + |
| Isobutanol | + | + |
| Isopropanol (2-Propanol) | + | + |
| Lactic acid | + | + |
| LB media | + | + |
| McCoy's 5A | + | + |
| MEM | + | + |
| Methanol | + | + |
| Methyl benzoate | | + |
| Methyl ethyl ketone | | + |
| Methyl propyl ketone | | + |
| Mineral oil (Engine oil) | | + |
| Monochloroacetic acid | | + |
| Nitric acid, 10% | | + |
| Octoxinol 9 (TRITON™ X-100) | + | + |
| Oxalic acid | + | + |
| PBS buffer | + | + |

| Reagent | seripettor® | seripettor® pro |
|--|-------------|--------------------|
| Perchloric acid | | + |
| Phenol | | + |
| Phosphoric acid, 85% | | + |
| Piperidine | | + |
| Polysorbate (TWEEN®) | + | + |
| Potassium chloride | + | + |
| Potassium dichromate | + | + |
| Potassium hydroxide | + | + |
| Potassium hydroxide in ethanol | + | + |
| Potassium permanganate | + | + |
| Propionic acid | + | + |
| Propylene glycol (Propanediol) | + | + |
| Pyridine | | + |
| Pyruvic acid | + | + |
| Ringer's solution | + | + |
| RPMI 1640 | + | + |
| Salicylaldehyde | | + |
| Salicylic acid | + | + |
| SDS (sodium dodecyl sulfate) | + | + |
| Silver acetate | + | + |
| Silver nitrate | + | + |
| Sodium acetate | + | + |
| Sodium chloride | + | + |
| Sodium dichromate | + | + |
| Sodium fluoride | + | + |
| Sodium hydroxide, 30% | + | + |
| Sodium hypochlorite 20% (active chlorine approx. 10%) | | + |
| Sulfuric acid, 10% | + | + |
| Tartaric acid | | + |
| Tris-buffered saline w. Tween20 | + | + |
| TE buffer | + | + |
| TRIS buffer | + | + |
| Urea | + | + |
| Zinc chloride, 10% | + | + |
| Zinc sulfate, 10% | + | + |

Note: seripettor® and seripettor® pro are not suitable for HF.

The above recommendations reflect testing completed prior to publication. Always follow instructions in the operating manual of the instrument as well as the reagent manufacturer's specifications. In addition to these chemicals, a variety of organic and inorganic saline solutions (e.g., biological buffers), biological detergents and media for cell culture can be dispensed. Should you require information on chemicals not listed, please feel free to contact BRAND. Status as of: 0420/9

BRAND®, seripettor®, BRAND. For lab. For life.® as well as the BRAND figurative mark are registered trademarks or trademarks of BRAND GMBH + CO KG, Germany. All other trademarks mentioned or depicted here are the property of the respective owners. Our technical literature is intended to inform and advise our customers. However, the validity of general empirical values, and of results obtained under test conditions, for specific applications depends on many factors beyond our control. Please appreciate, therefore, that no claims can be derived from our advice. The user is responsible for checking the appropriateness of the product for any particular application.

California Residents: For more information concerning California Proposition 65, please refer to www.brand.de/calprop65
Subject to technical modification without notice. Errors excepted.

BRAND GMBH + CO KG · P.O. Box 11 55 · 97861 Wertheim · Germany
Phone: +49 9342 808-0 · Fax: +49 9342 808-98000 · E-Mail: info@brand.de · Internet: www.brand.de

