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➔ DANGER indicates a hazardous situation which, if not avoided, will result in death or serious injury.

+ WARNING indicates a hazardous situation which, if not avoided, could result in death or serious injury.

• CAUTION indicates a hazardous situation which, if not avoided, could result in minor or moderate injury.

NOTICE is used to address practices not related to personal injury.

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After sales service:
Contact your local dealer or call +49 9342 808-5500.

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Trademark index:

VACUU•LAN® (US-Reg.No 3,704,401), VACUU•BUS®, VACUU•CONTROL®, Peltronic®, VARIO® (US-Reg.No 3,833,788), VACUUBRAND® (US-Reg.No 3,733,388) and also the shown company logos are registered trademarks of VACUUBRAND GMBH + CO KG in Germany and/or other countries.

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Safety information!

DANGER

- ➔ Ensure that the pump cannot be operated accidentally. Never operate the pump if covers or other parts of the pump are disassembled.
- ➔ Before starting work, disconnect the electrical power cord.
- ➔ **Note: The pump might be contaminated with the process chemicals that have been pumped during operation. Ensure that the pump is decontaminated before maintenance. Avoid the release of pollutants.**

WARNING

- ☞ **Read the pump's manual carefully before installing or operating the equipment. Observe the instructions contained in the pump's manual.**
- ☞ Ensure that the individual components are only connected, combined and operated according to their design and as indicated in the instructions for use. Use only **original manufacturer's spare parts and accessories**. Otherwise the safety and performance of the equipment, as well as the electromagnetic compatibility of the equipment might be reduced.
The CE mark or the cTÜVus mark may be voided if not using original manufacturer's spare parts.
- ☞ Never use defective or damaged parts. Replace defective or damaged parts in any case.
- ☞ Take adequate precautions to protect people from the effects of dangerous substances that may have contaminated the pump and may be released upon disassembly. Ensure that the maintenance technician is familiar with the safety procedures which relate to the products processed by the pumping system.
Use appropriate protective clothing, safety goggles and protective gloves.
- ☞ Vent the pump and isolate it from the vacuum system before you start maintenance. Allow sufficient cooling of the pump before starting work.
- ☞ **Attention:** Comply with applicable regulations when disposing of chemicals, condensates or oils. Take into consideration that those substances may be contaminated

NOTICE

This manual is intended for trained service people only. Ensure that work is done only by a suitably trained and supervised technician. Comply with local and national safety requirements.

Before starting work check that the required parts are available and of the correct type. Check that the required tools are available.

- Check the operating sequence at first mentally on feasibility, safety requirements and consequences on safety and function of the equipment.
- Check the pump's operability and safety after work.
- In case of problems occurring, return the equipment to the factory for inspection or repair, if necessary.

In order to comply with laws (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments can only be returned when certain procedures (see section **"Notes on return to the factory"** in the pump's instructions for use) are followed.

Scrapping and waste disposal:

Dispose of the equipment and any components removed from it safely in accordance with all local and national safety and environmental requirements. Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from your processes.

- You may authorize us to dispose of the equipment **at your expense**.

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Notes on using separators

NOTICE

Check the liquid level in the catchpots during operation. Check the liquid level in the catchpots regularly. Do not allow the catchpots to overflow.

Separator at the inlet:

Drain catchpot in good time, never allow liquid to reach the middle inlet pipe.

Oil mist separator:

Drain catchpot in good time, never allow liquid level to reach the filter cartridge.

Oil mist filter	Maximum permissible amount of liquid in the separator catchpot
Oil mist filter RE/RZ 2 / 2.5 / 5 / 6 (20698003)	max. 180 ml (till bottom edge of filter cartridge)
Oil mist filter RE/RZ 8 / 16 (20698008)	max. 250 ml (till bottom edge of filter cartridge)
Oil mist filter RE/RZ 9 / 16 (20698017)	max. 195 ml (till bottom edge of filter cartridge)

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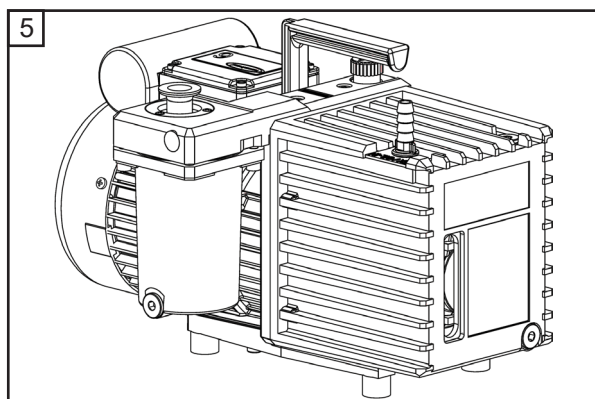
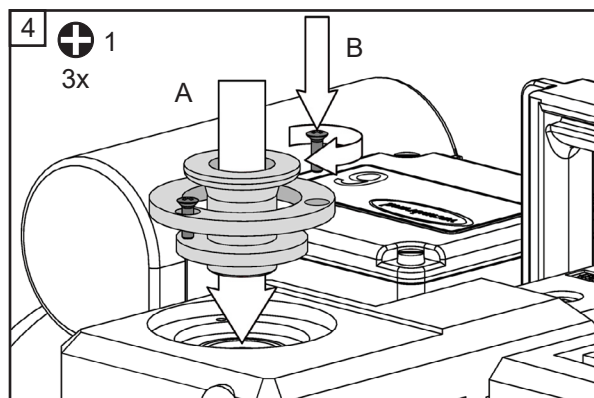
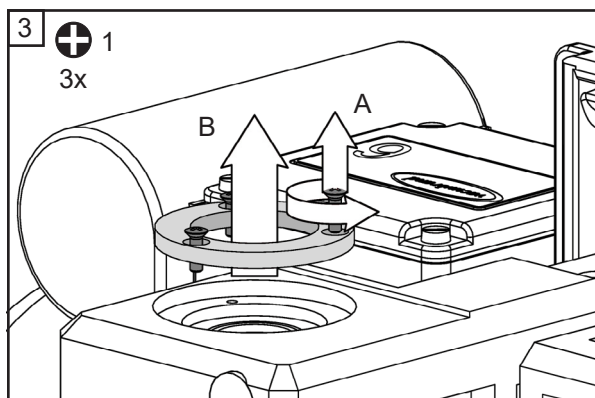
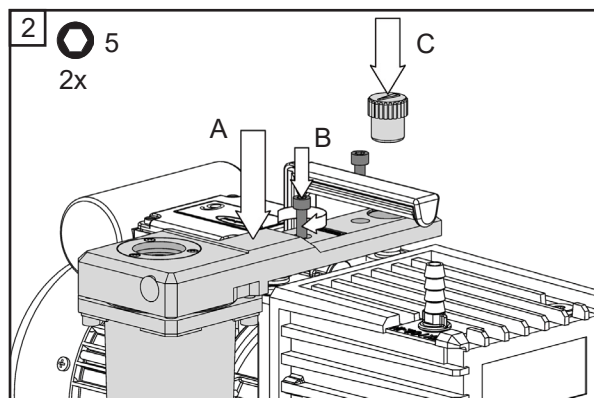
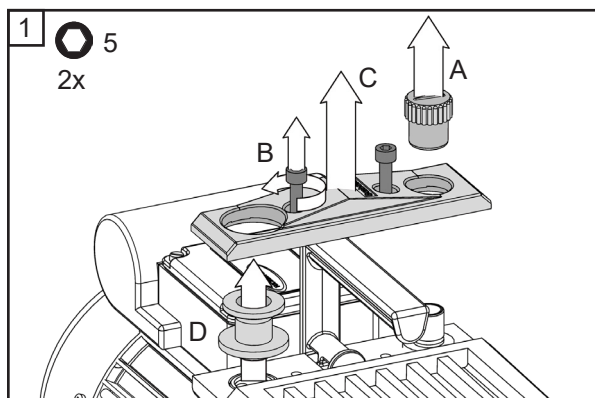
 Check that seals are correctly positioned and in perfect condition.

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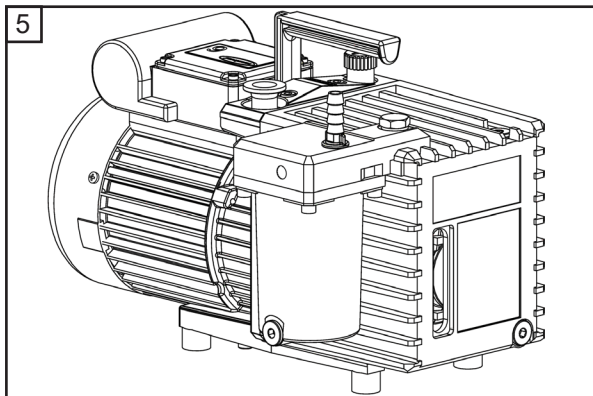
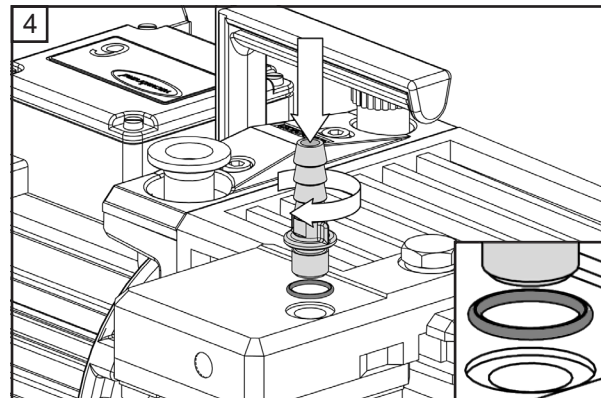
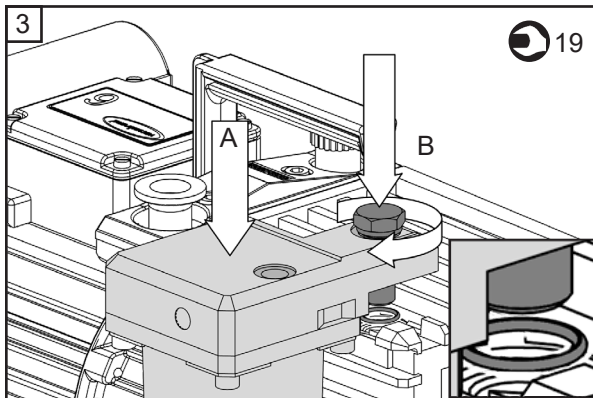
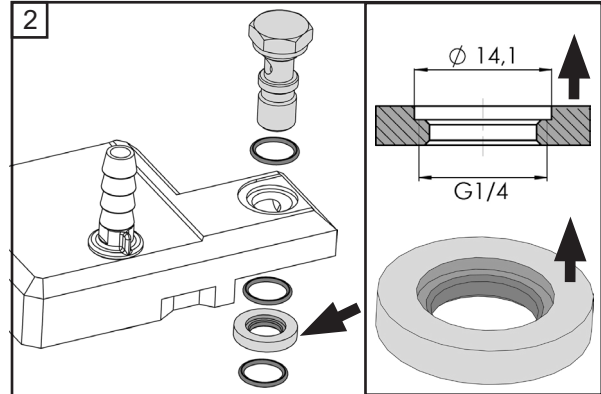
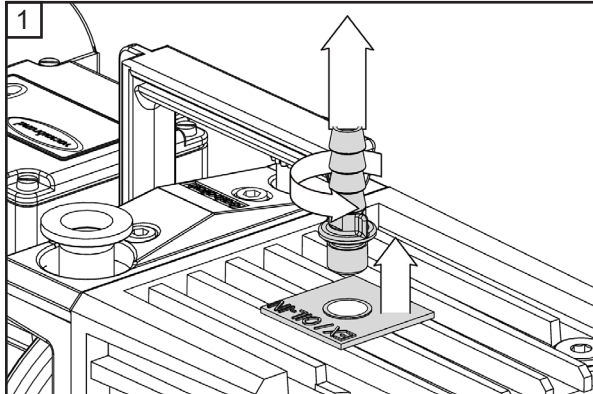
 **Note:** Aluminum thread. Aluminum parts can easily be damaged. Install screw with care and do not overtighten.





Wechsel des Filterelements im Auspufffilter / Replacing the filter element in the oil mist filter RE / RZ 2 / 2.5 / 5 / 6 (20698004)

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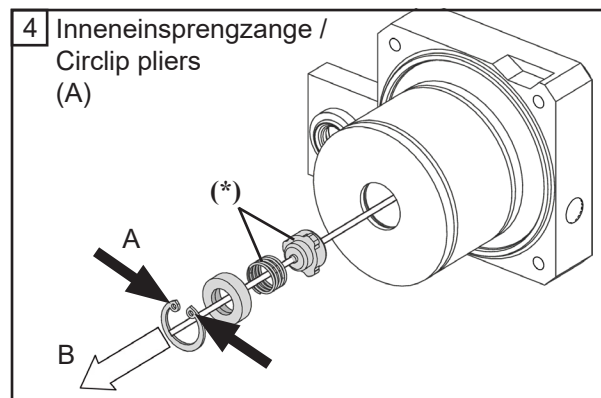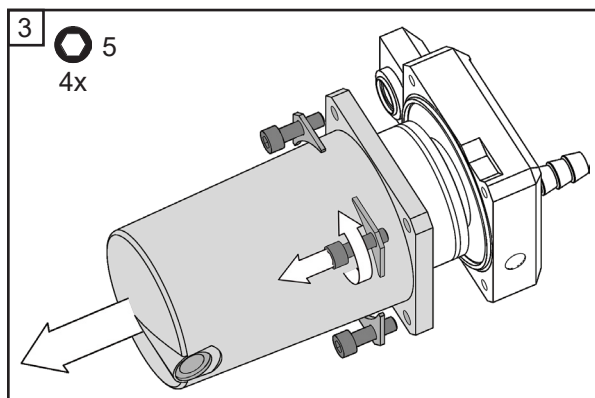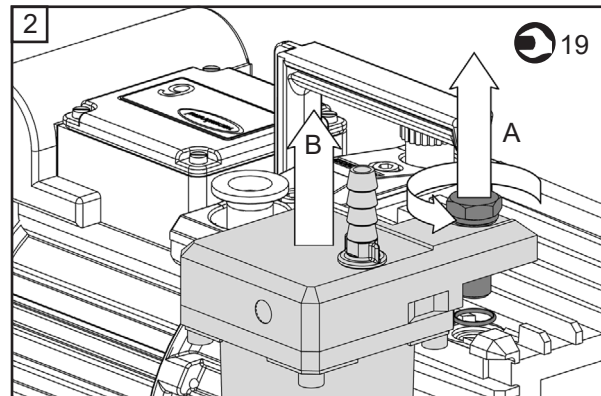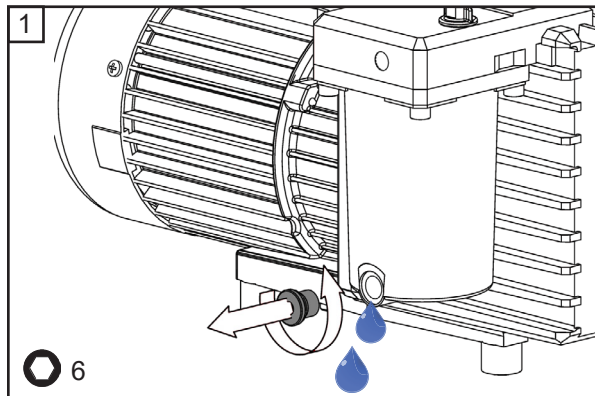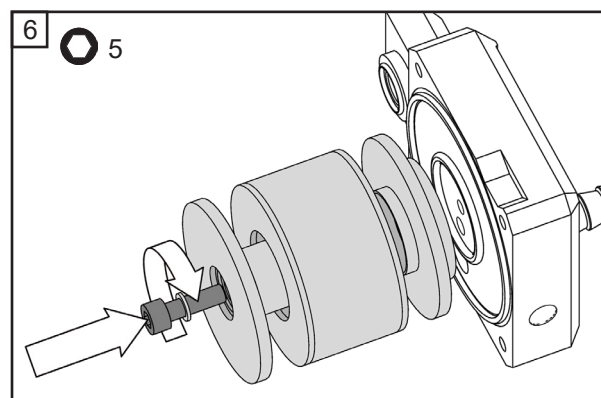
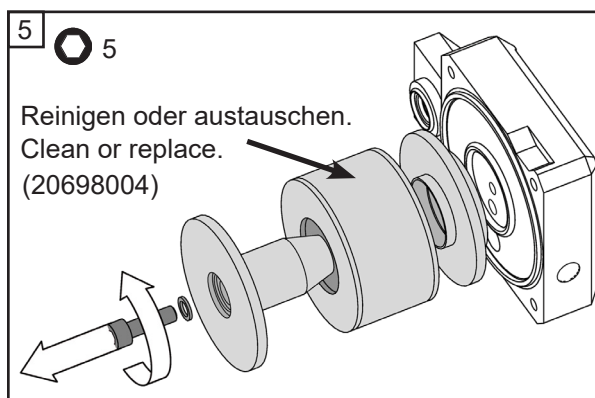
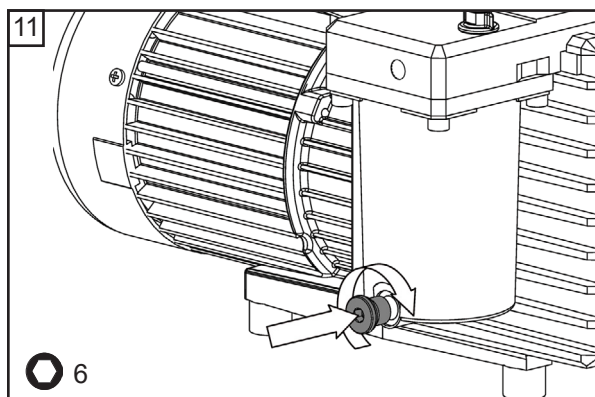
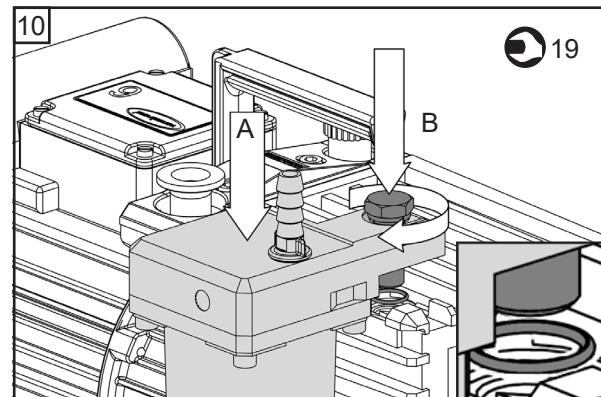
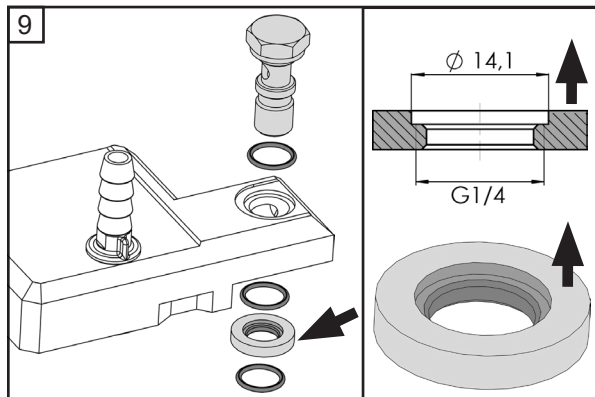
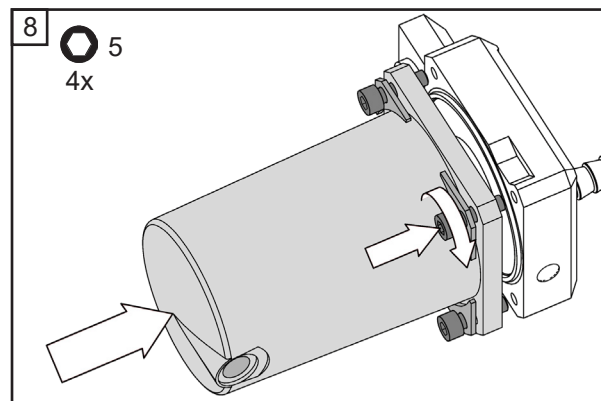
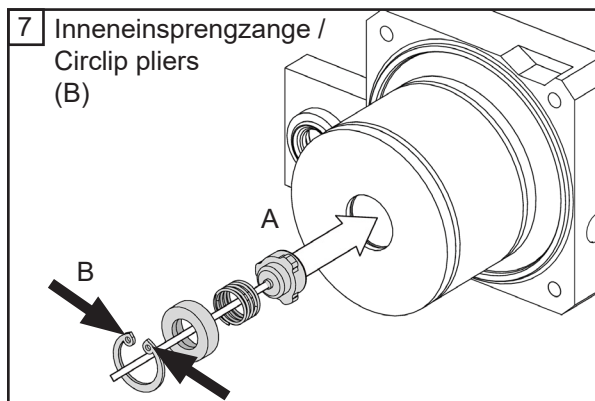
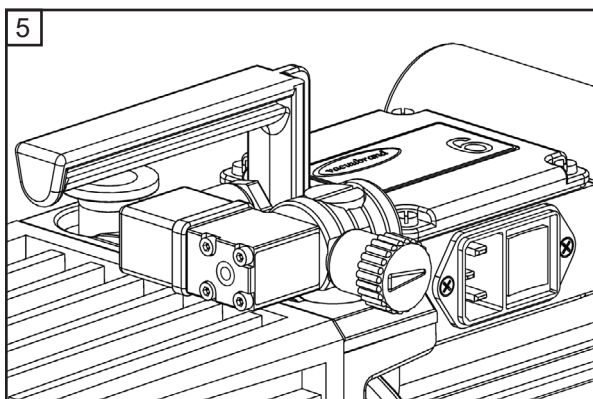
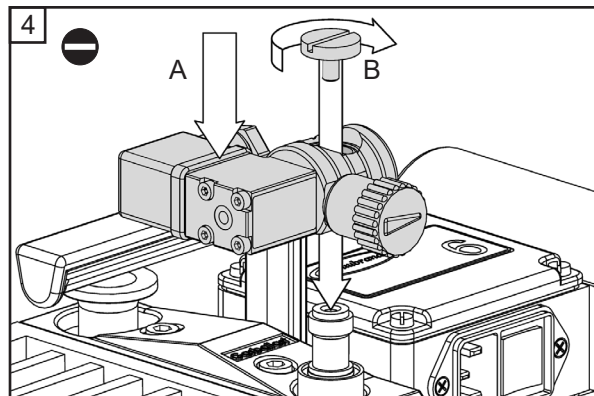
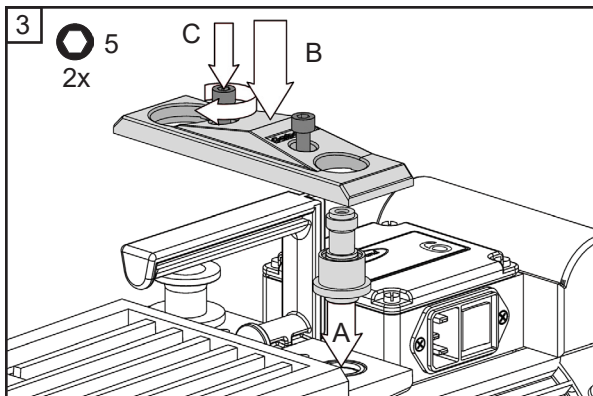
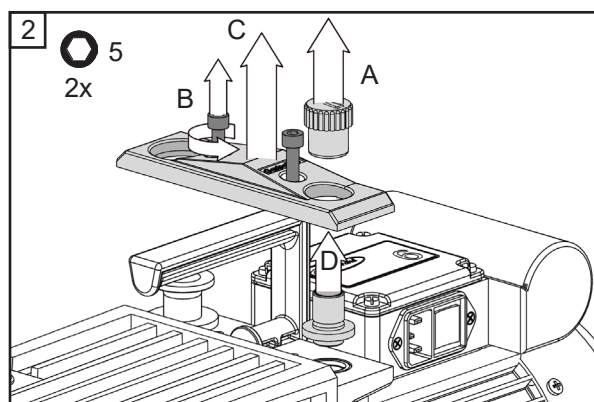
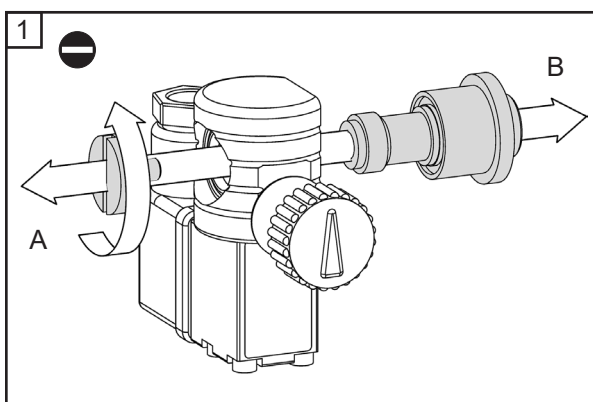
☞ Check that the O-ring (3 A) is correctly positioned and in perfect condition.

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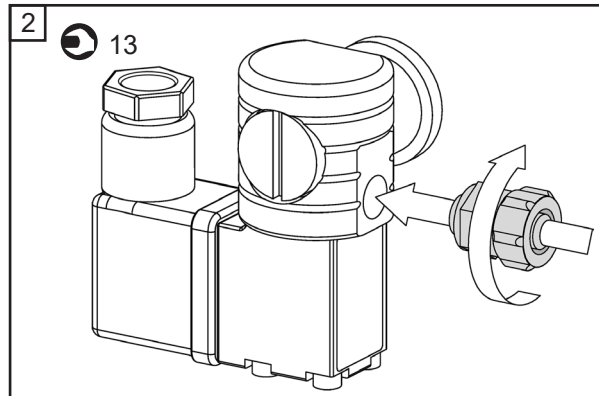
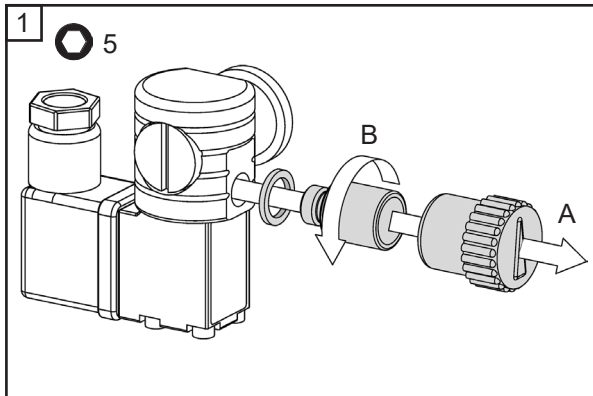
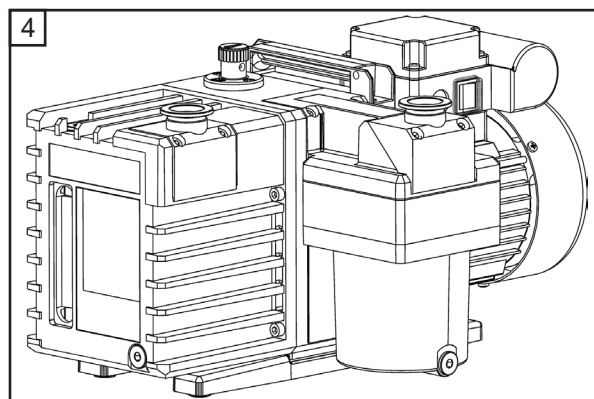
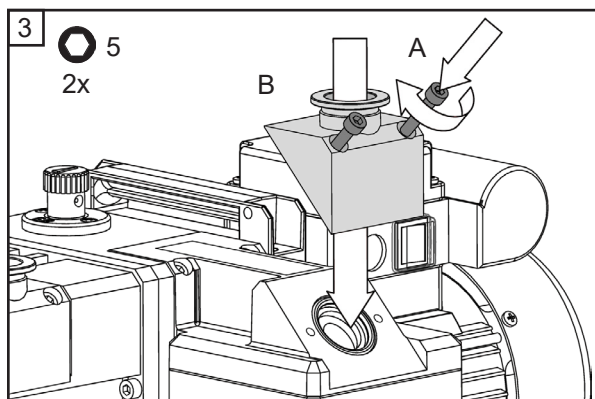
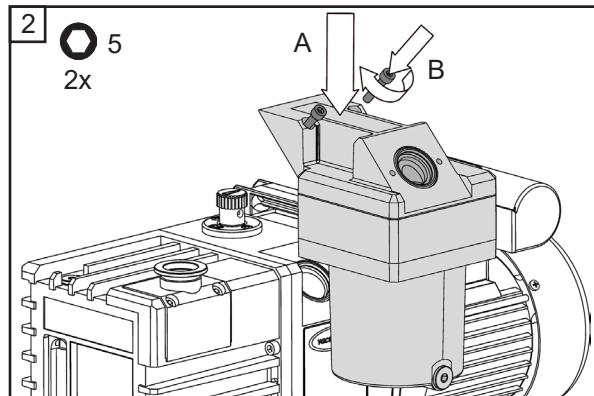
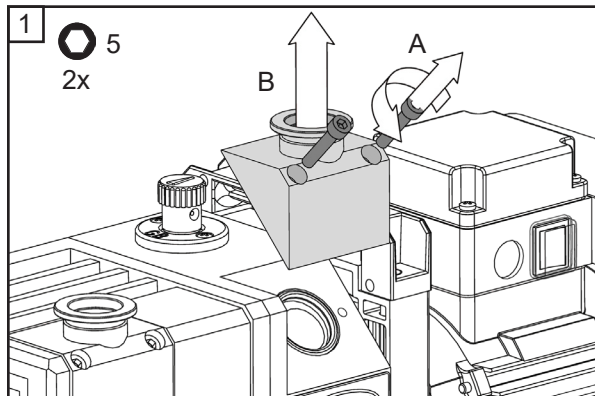
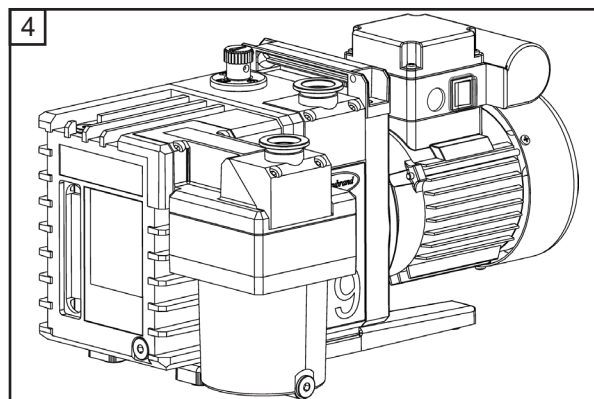
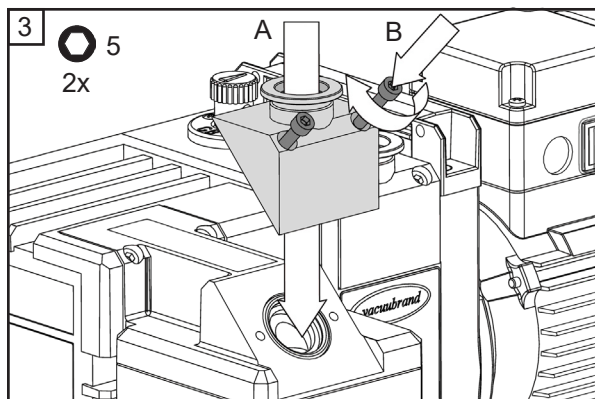
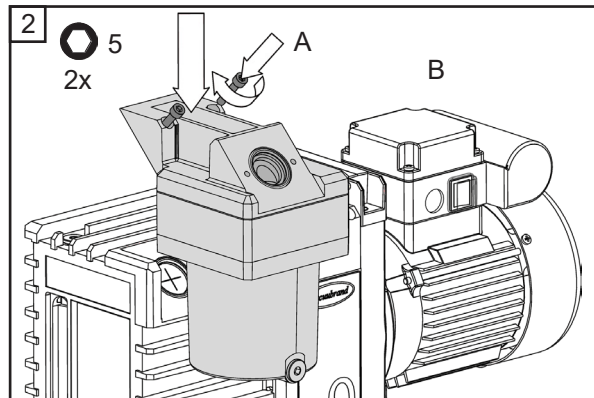
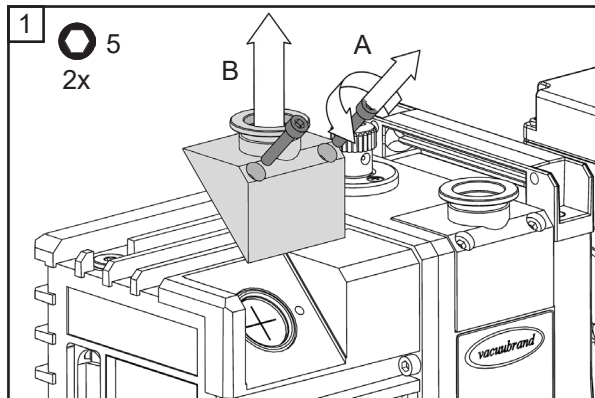
☞ Check that seals are correctly positioned and in perfect condition.







Wechsel des Filterelements im Auspufffilter (20698017) / Replacing the filter element in the oil mist filter (20698017) RE / RZ 9 / 16 (20667085)

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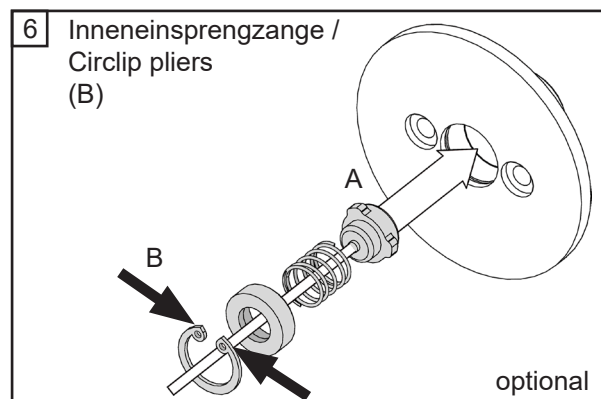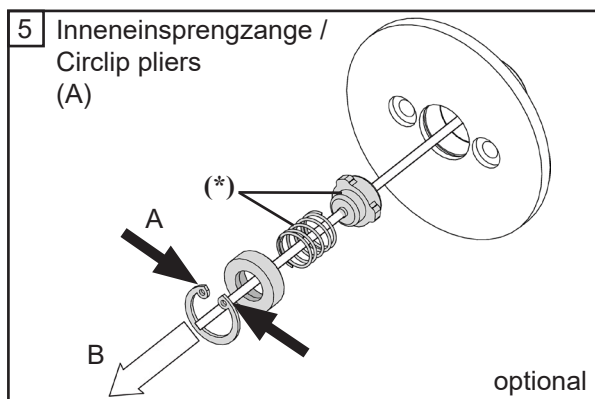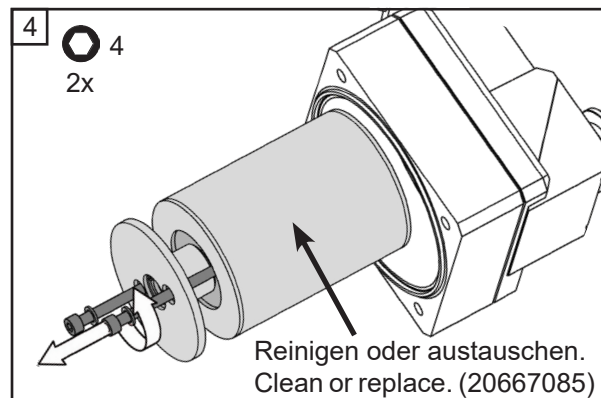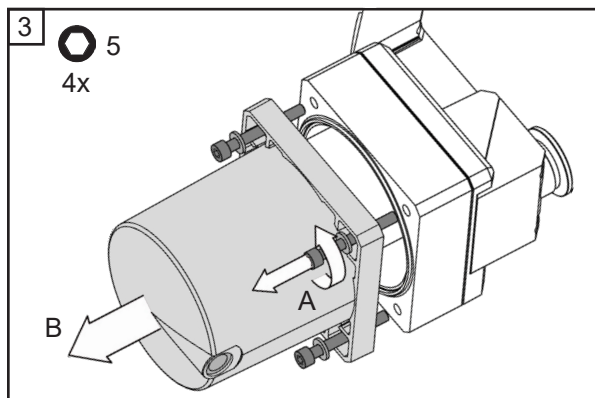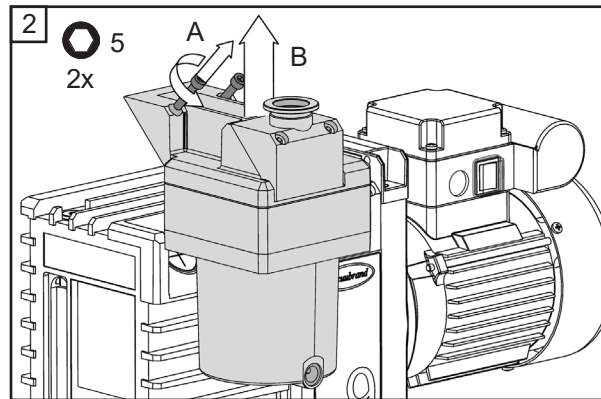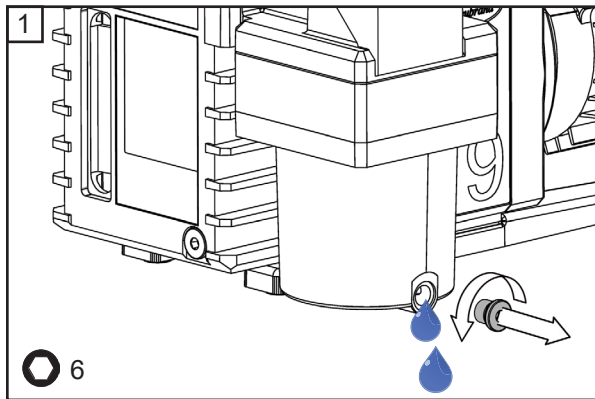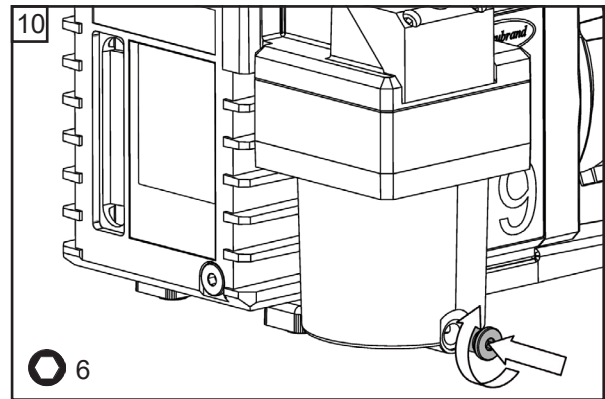
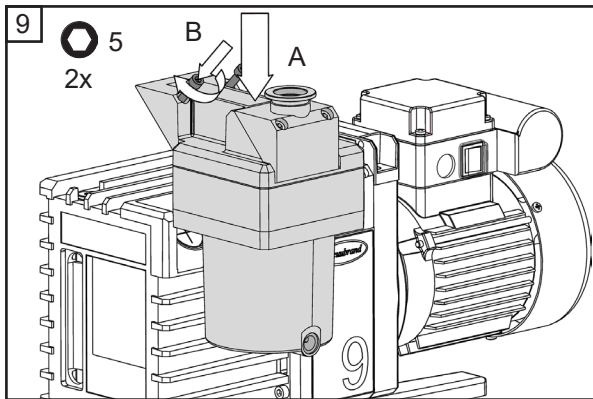
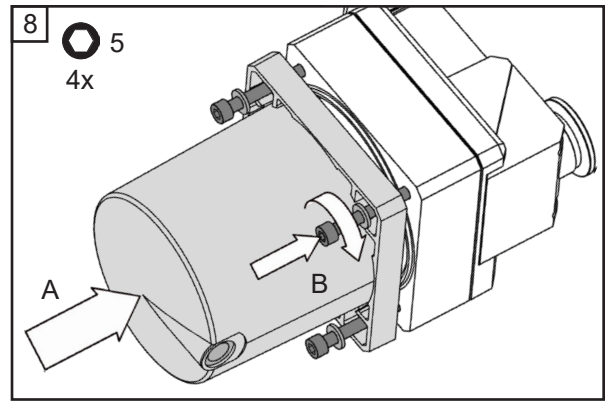
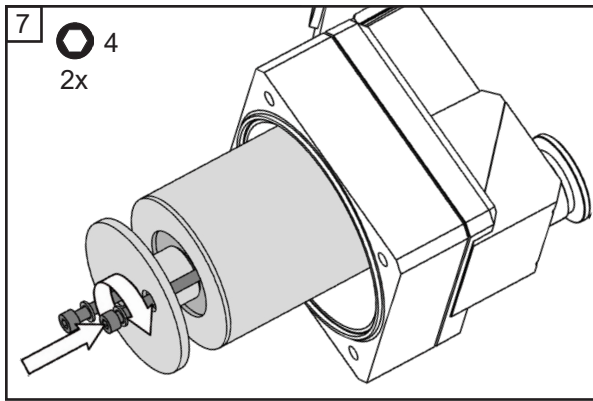
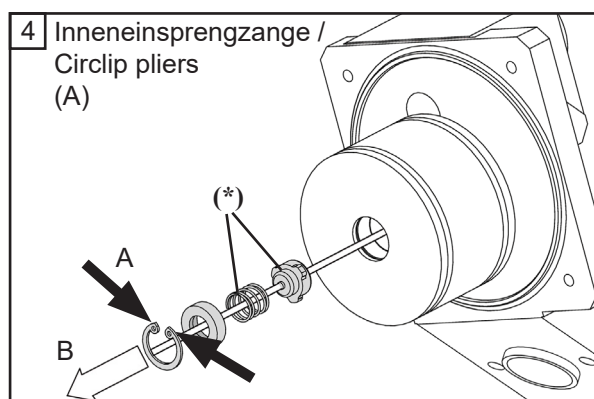
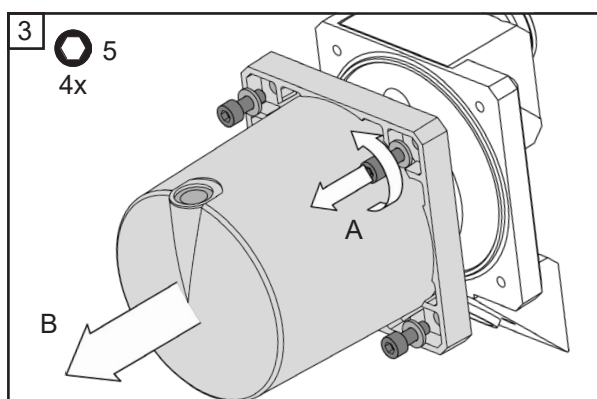
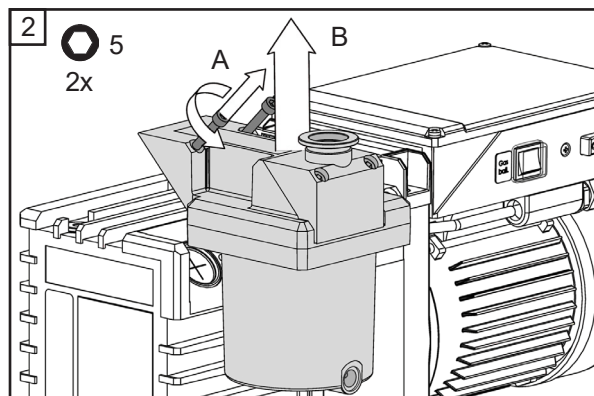
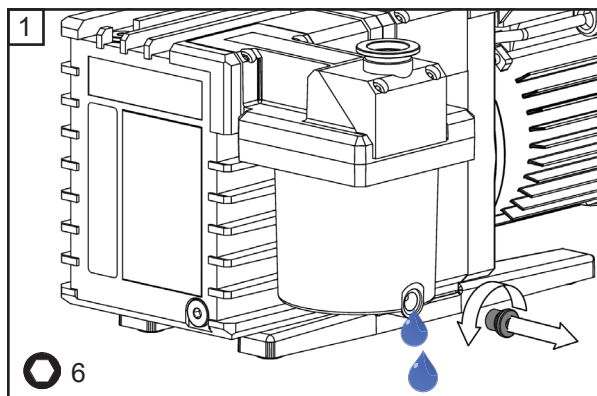
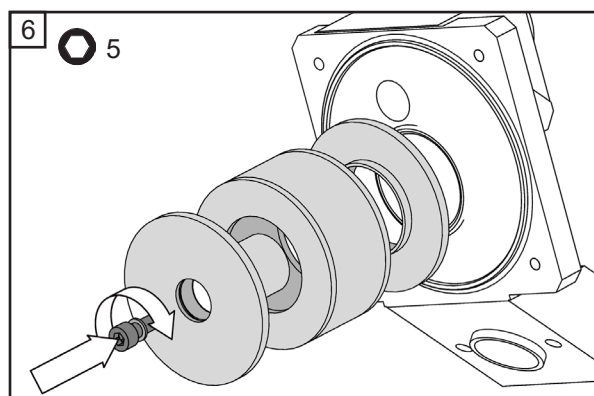
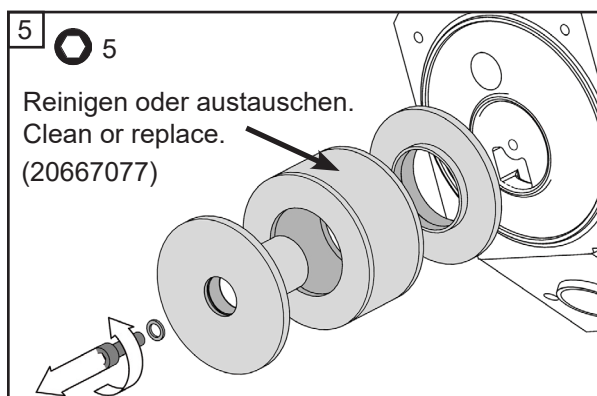
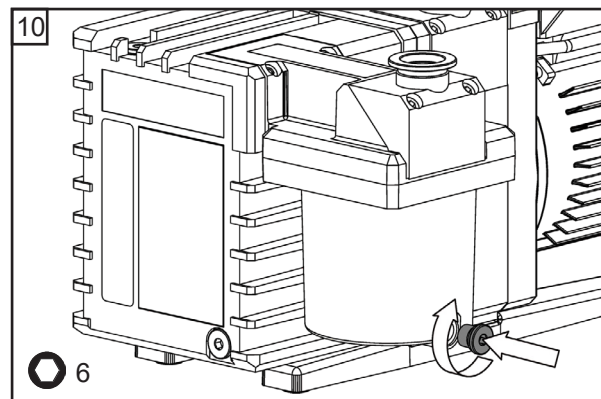
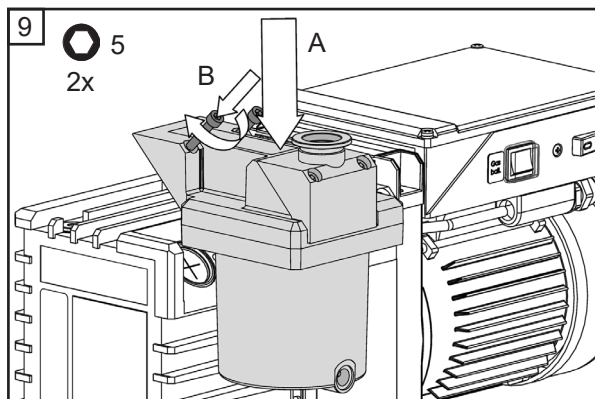
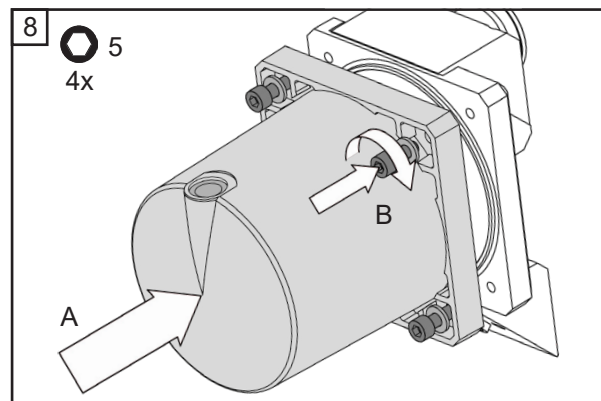
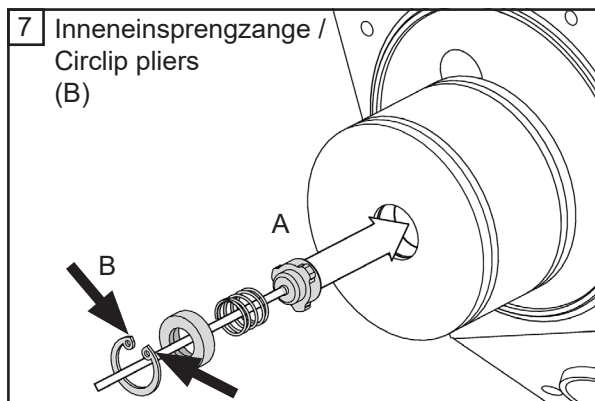
☞ Check that valve spring (*) and valve cap (*) are in perfect condition.



Hauptstromölfilter / Full-flow oil filter RE / RZ 8 / 9 / 16 (20698010)

Achtung: Hauptstromölfilter nur an Pumpen bis Baujahr 06/2016 montierbar.
Attention: Full-flow oil filter fits only for pumps until year of manufacture 06/2016.

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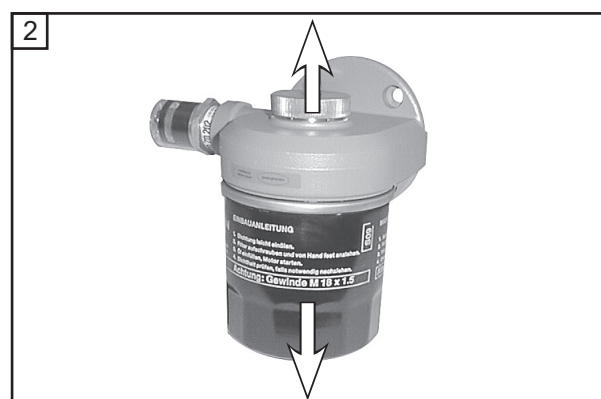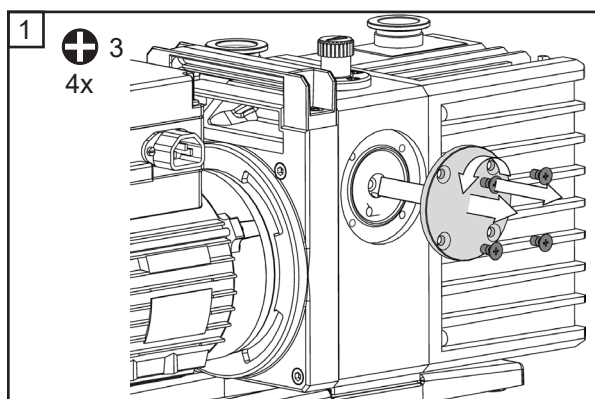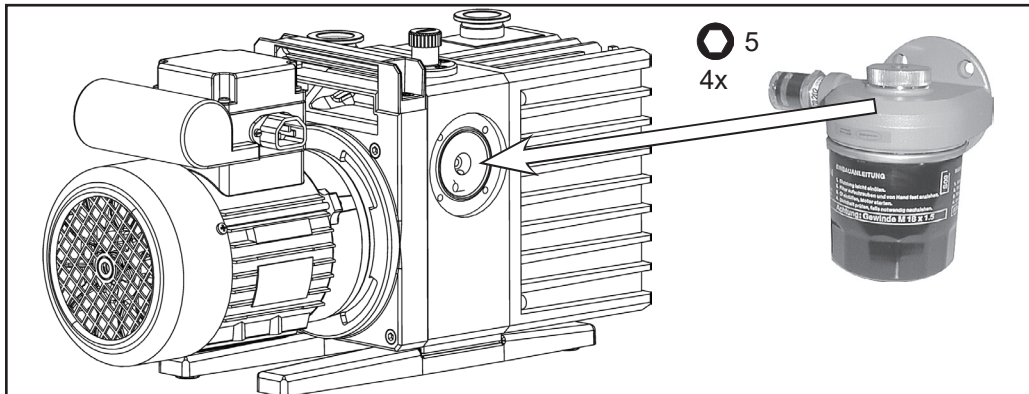
- ➔ Fix housing of full flow oil filter with four Allen screws to pump housing.
- ➔ Pour new pump oil into filter element.
- ➔ Coat seals of filter with oil and install filter. Tighten until hand-tight.
- ➔ Switch on pump and check filter for leaks. Tighten filter if necessary.
- ➔ Check oil level. Top up if necessary.

Wechsel des Hauptstromölfilters / Replacing the full-flow oil filter **RE / RZ 8 / 9 / 16 (20698010)**

Achtung: Hauptstromölfilter nur an Pumpen bis Baujahr 06/2016 montierbar.
Attention: Full-flow oil filter fits only for pumps until year of manufacture 06/2016.

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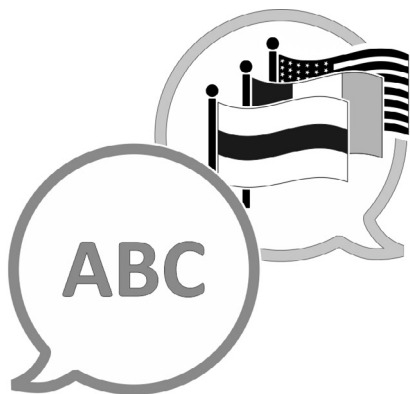
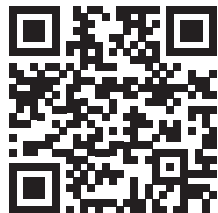
- ☞ Important: Oil might drip from pump - make sure surface beneath pump is suitable.
- ➔ Fix the lower part of the filter element by hand and unscrew banjo bolt (8 mm wide Allen key) at the top of the filter. Remove filter element.
- ☞ **Note:** Filter contains oil. Dispose of oil and filter in accordance with regulations.
- ➔ Coat seals of new filter with oil and install filter. Tighten until hand-tight.
- ➔ Switch on pump and check filter for leaks. Tighten filter if necessary.
- ➔ Check oil level. Top up if necessary.

[VACUUBRAND > Support > Manuals](#)

Manufacturer:

VACUUBRAND GMBH + CO KG
Alfred-Zippe-Str. 4
97877 Wertheim
GERMANY

Phone:

Head office +49 9342 808-0

Sales +49 9342 808-5550

Service +49 9342 808-5660

Fax: +49 9342 808-5555

Email: info@vacuubrand.com

Web: www.vacuubrand.com